NOVINKY A VYCHYTÁVKY V CHROMATOGRAFII THERMO SCIENTIFIC

Lukáš Plaček
Michaela Rajčanová

pragolab

ThermoFisher
SCIENTIFIC

Iontová Chromatografie

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

CHROMELEON 7.2
Simply Intelligent

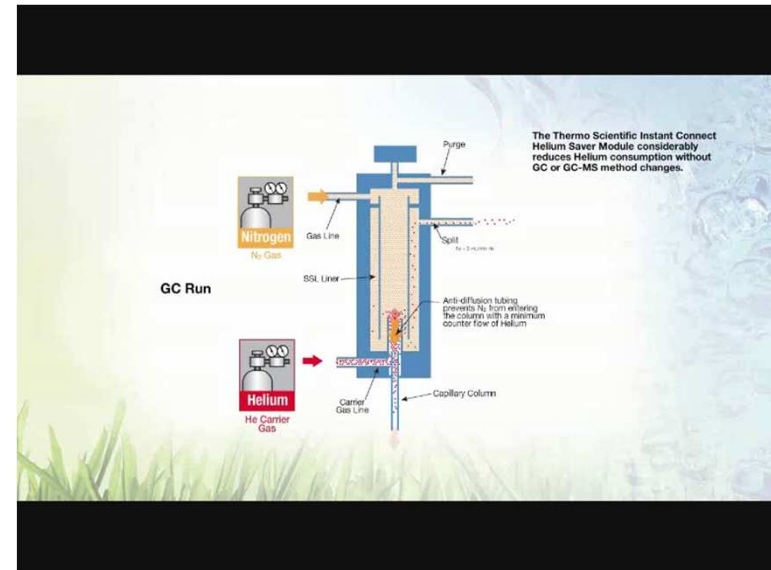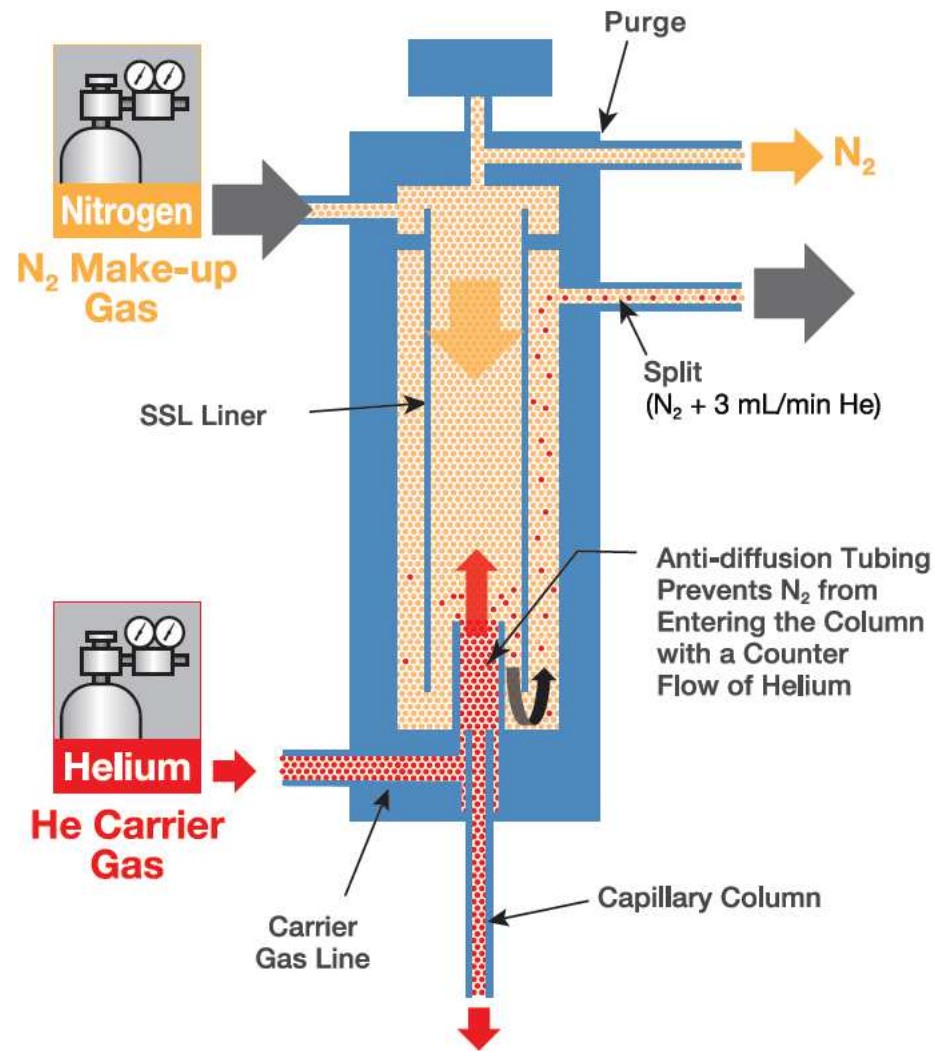GC: helium saver modul

Helium Saver Extends the Lifetime of the Helium Cylinder

3.5 years continuously used 24/7/365 for GC-MS analysis

5.4 years shutting helium off or to N₂ during weekends

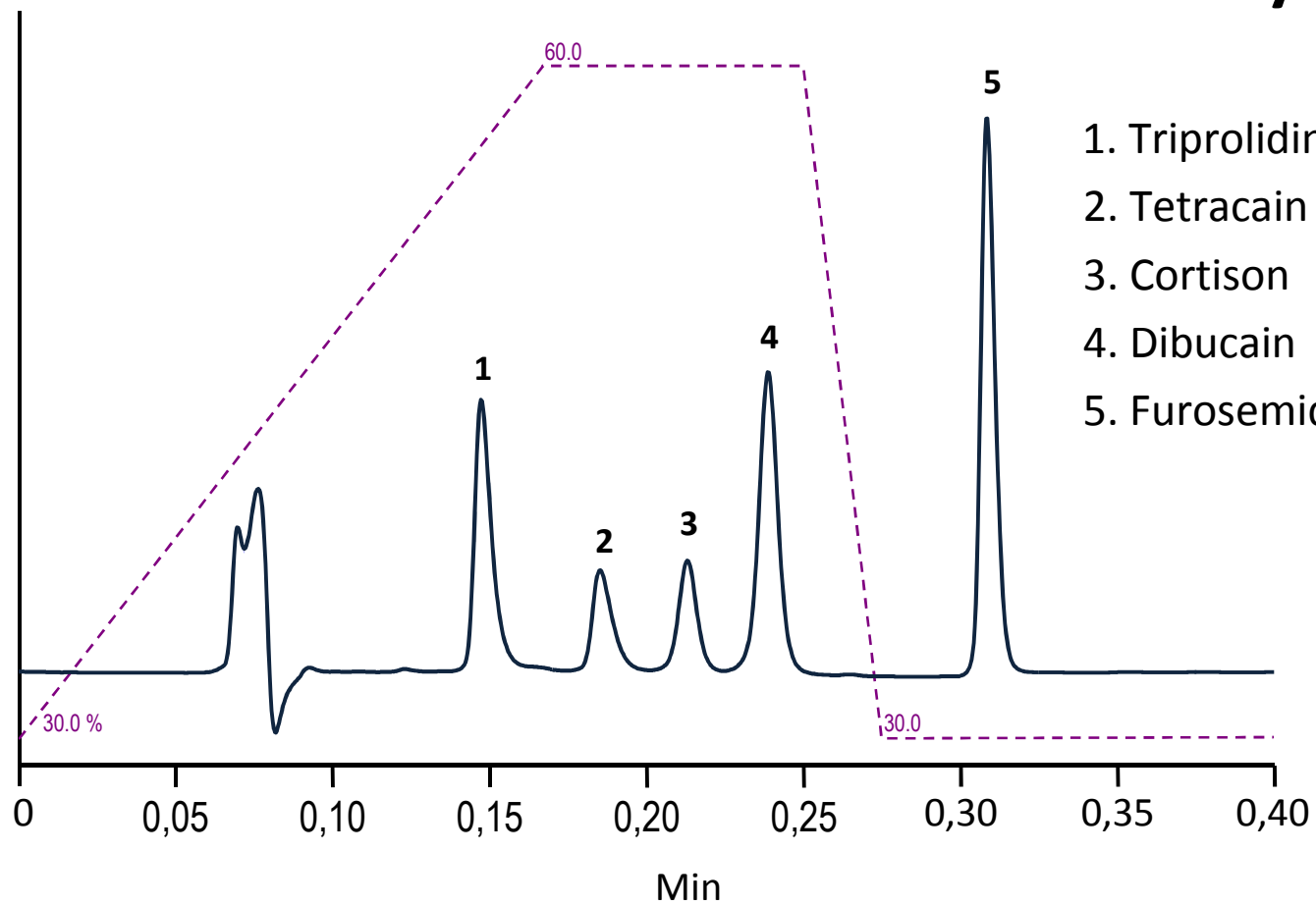
10.5 years shutting helium off or to N₂ overnight (8 hr day)

14.6 years shutting helium off or to N₂ on weekends and overnight

This could be the **last and only** helium cylinder that will be needed for the lifetime of the instrument

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
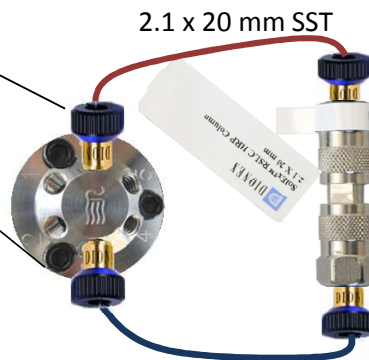
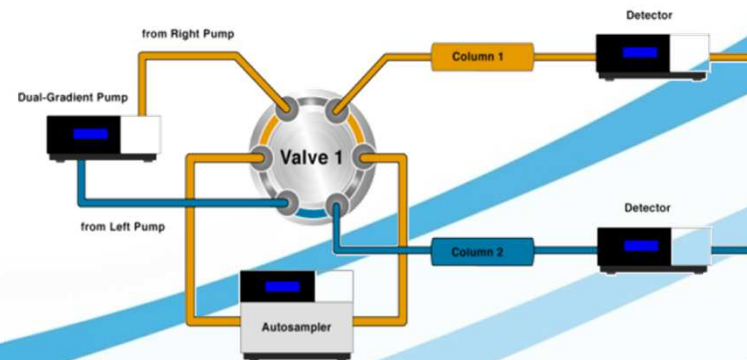
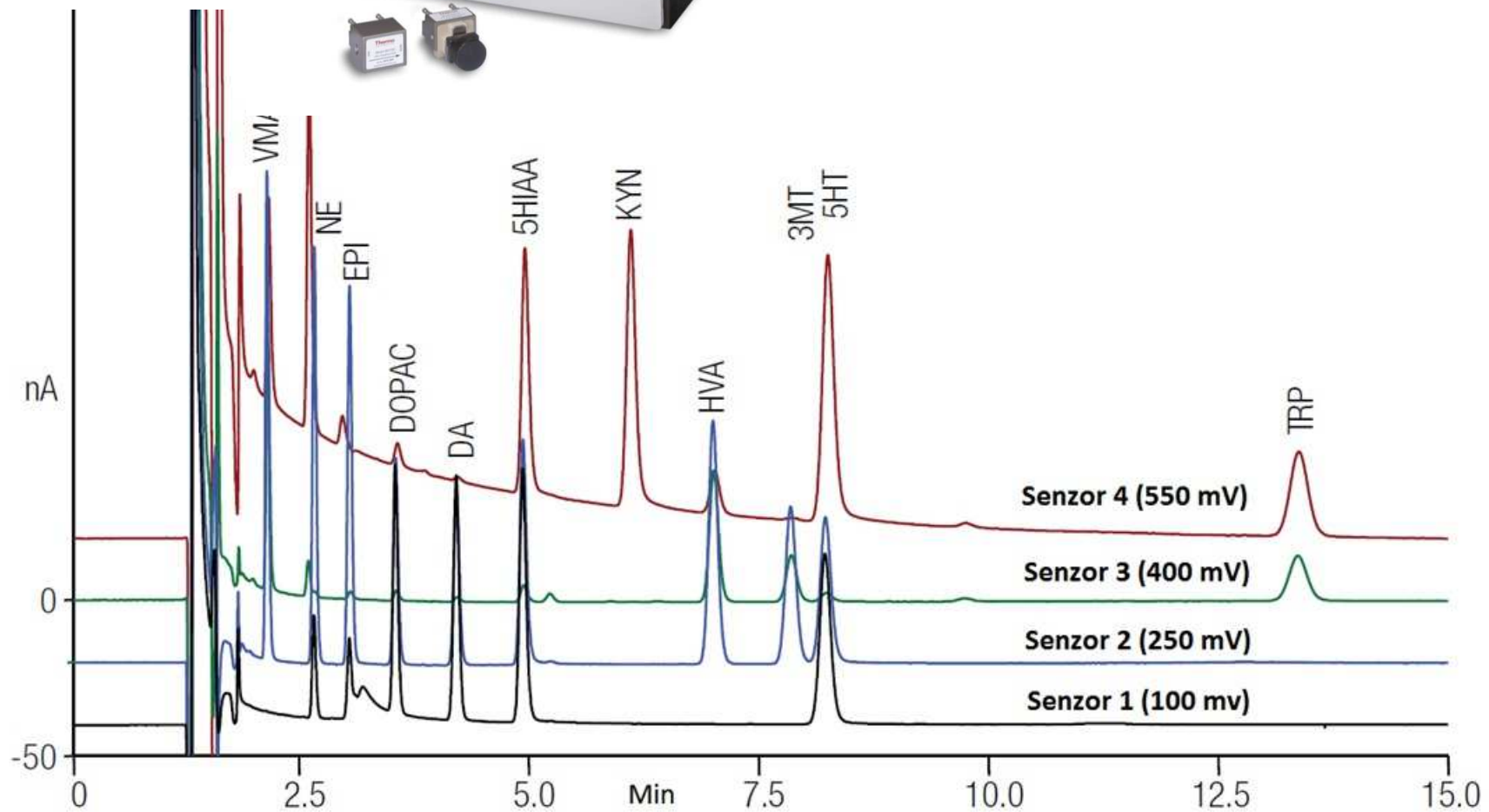
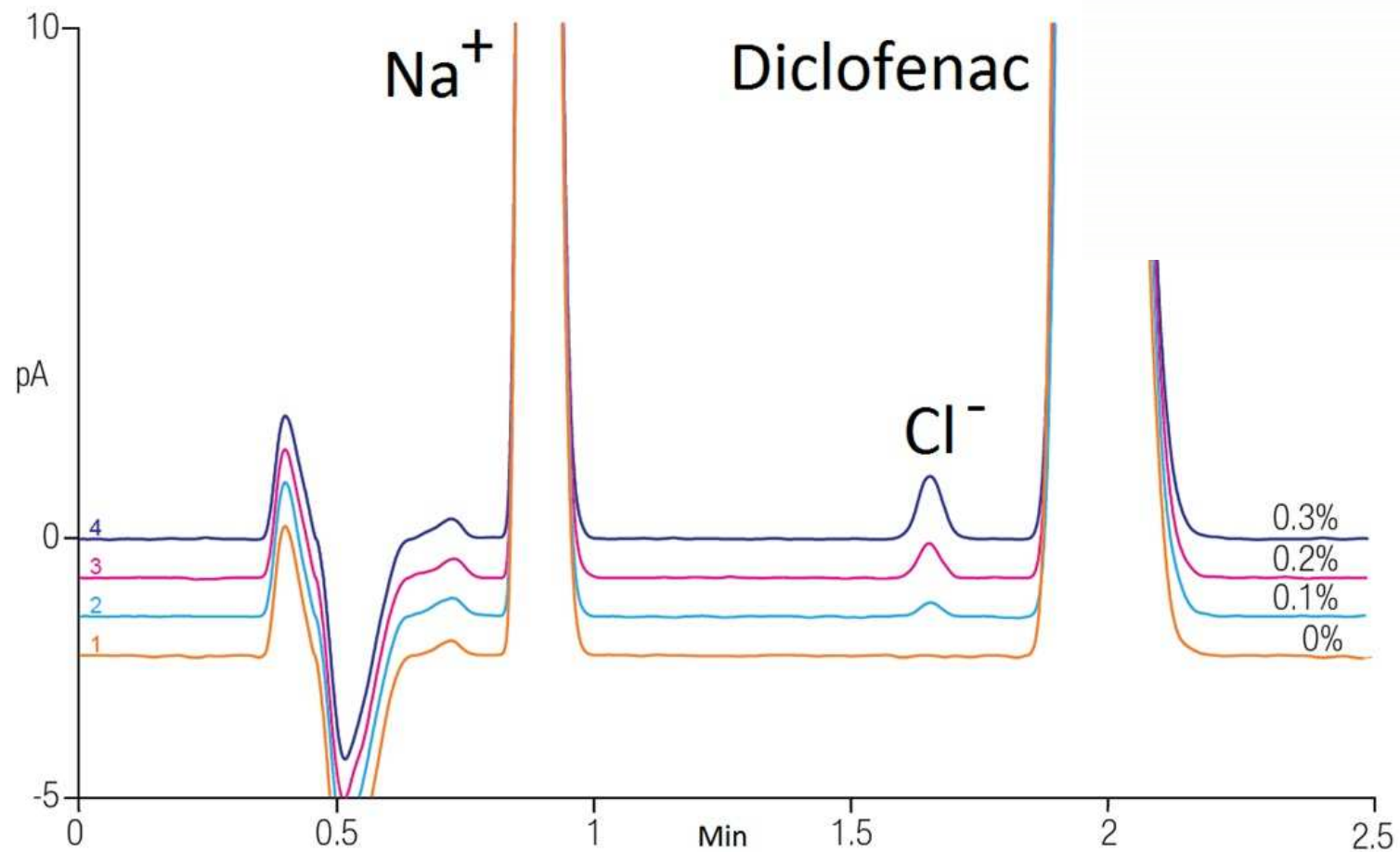
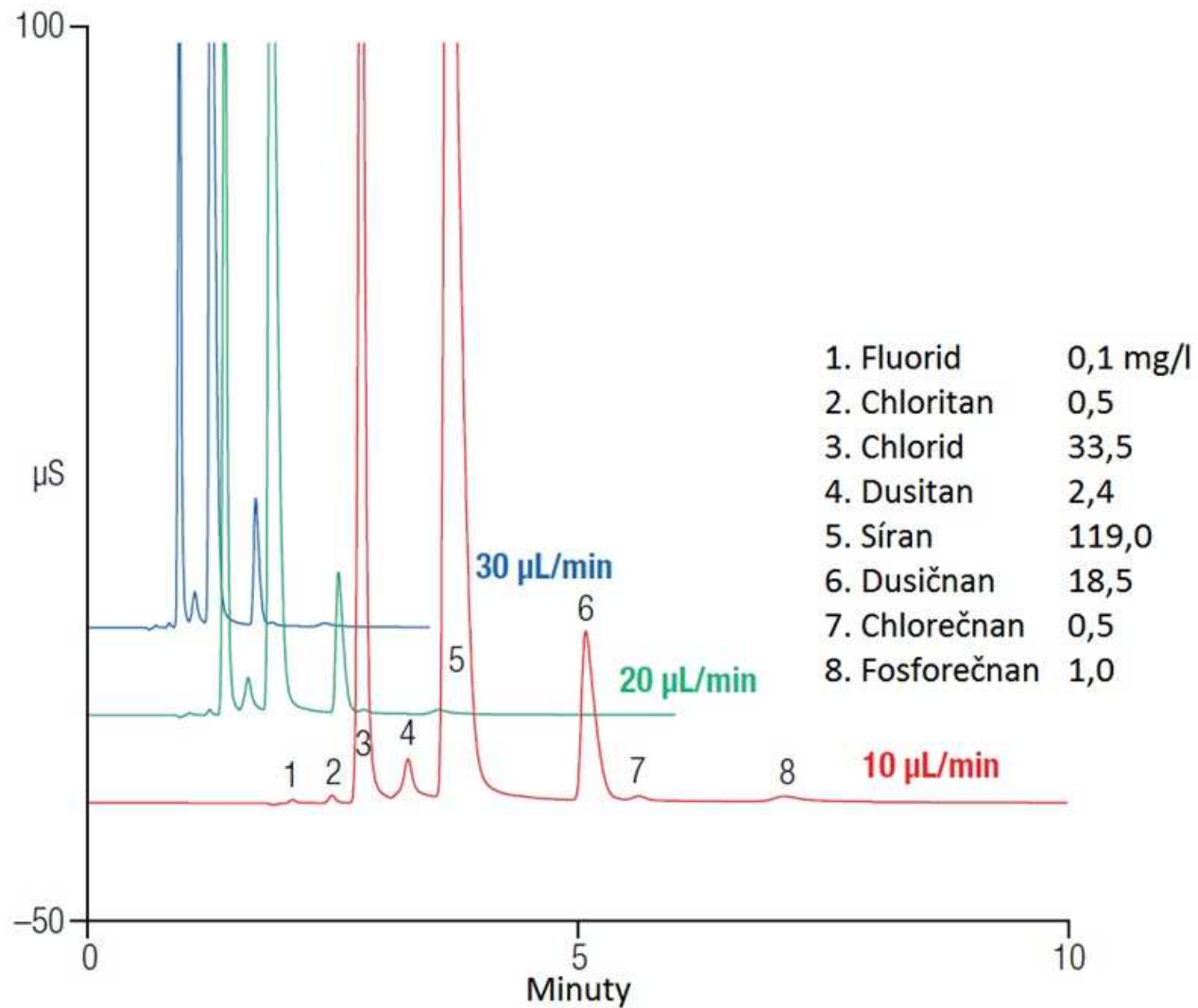
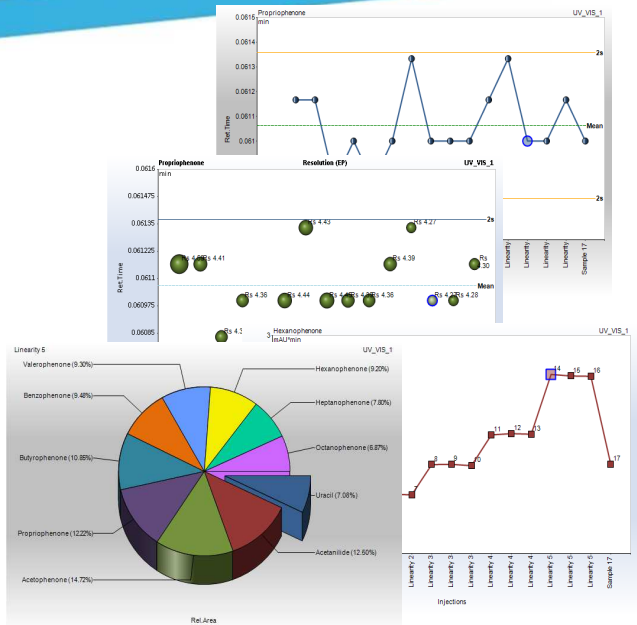
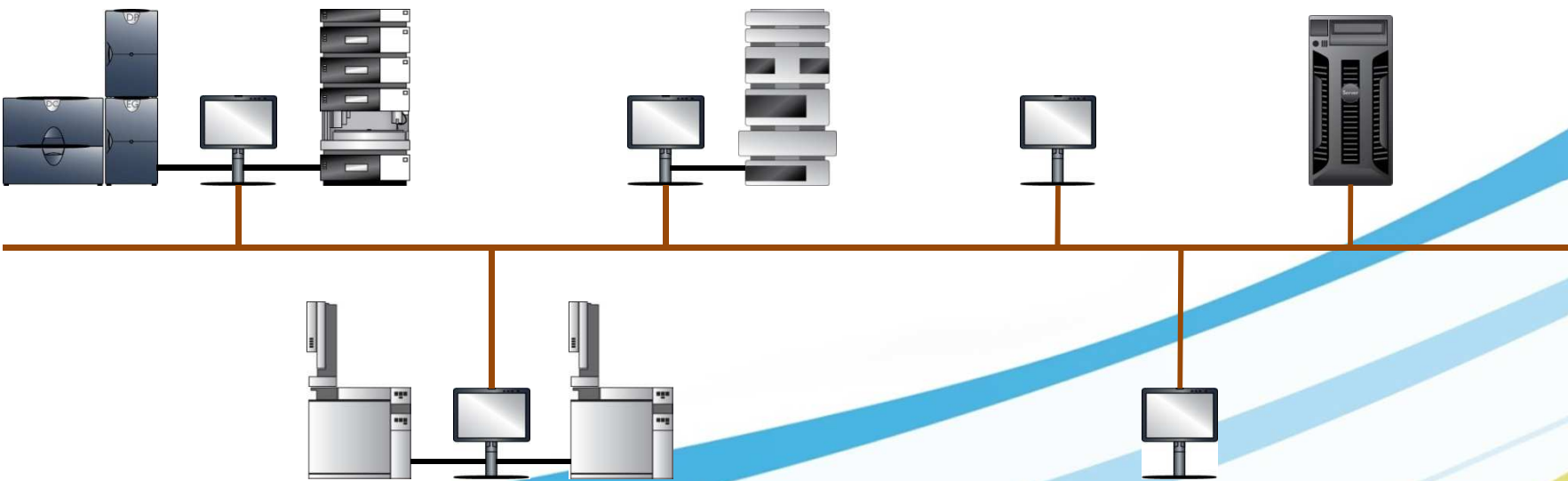
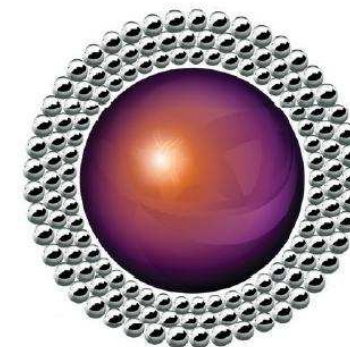
- ✓ 2,2 μm


Nano-polymer bead (SAX)

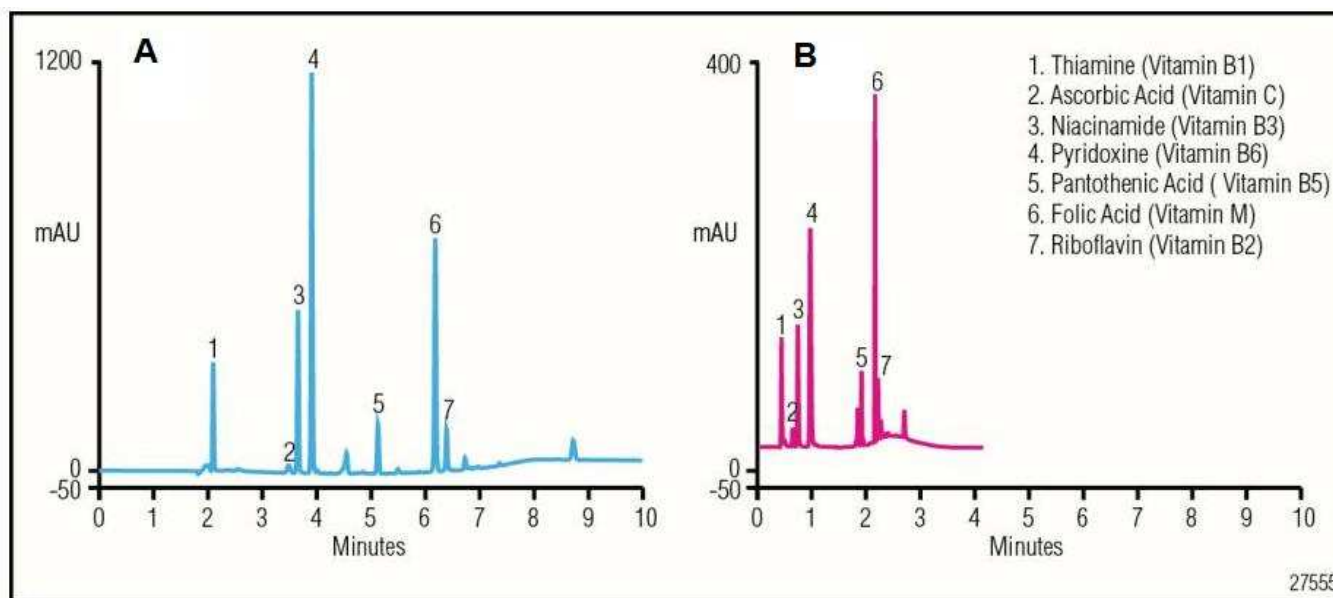Analysis of seven water-soluble vitamins under A) conventional LC and B) UHPLC conditions on Acclaim PA2, 5 μm , 250 \times 4.6 mm and Acclaim RSLC PA2, 2.2 μm , 100 \times 2.1 mm columns, respectively.

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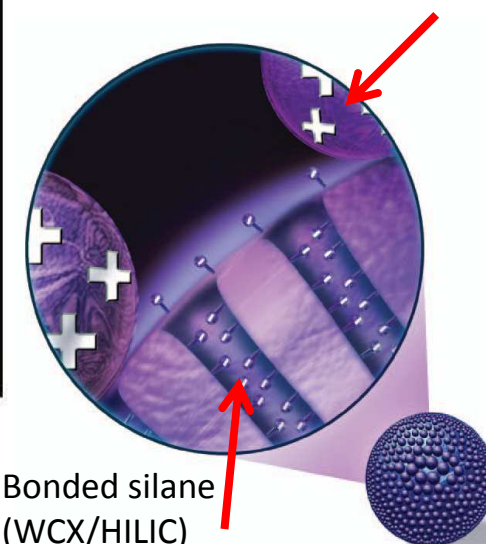
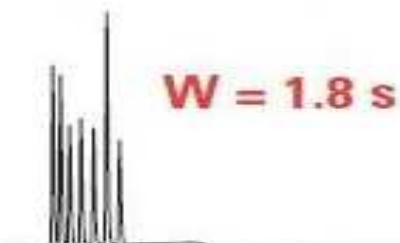
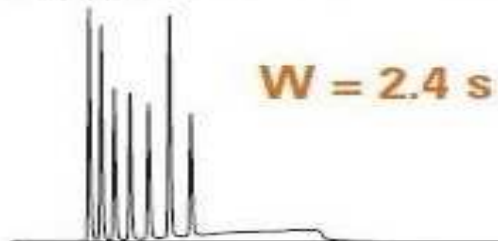
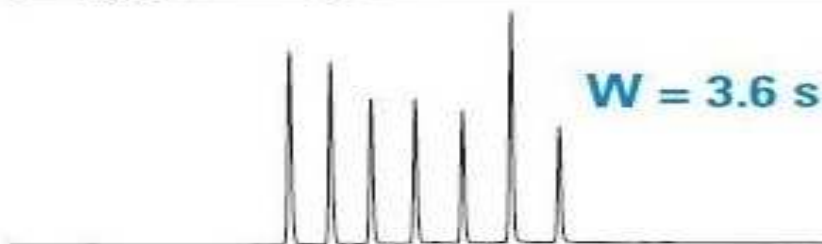
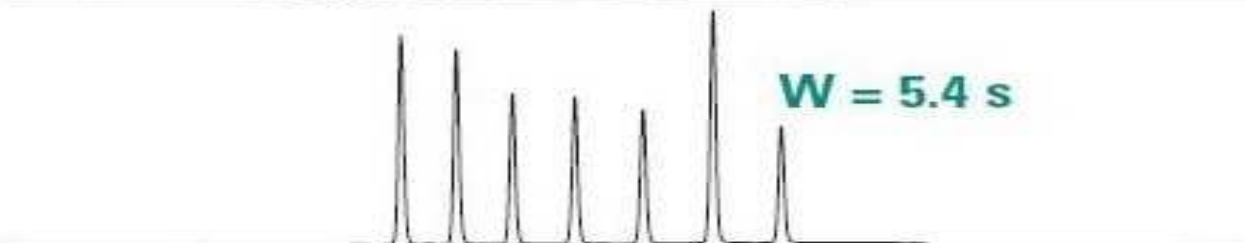
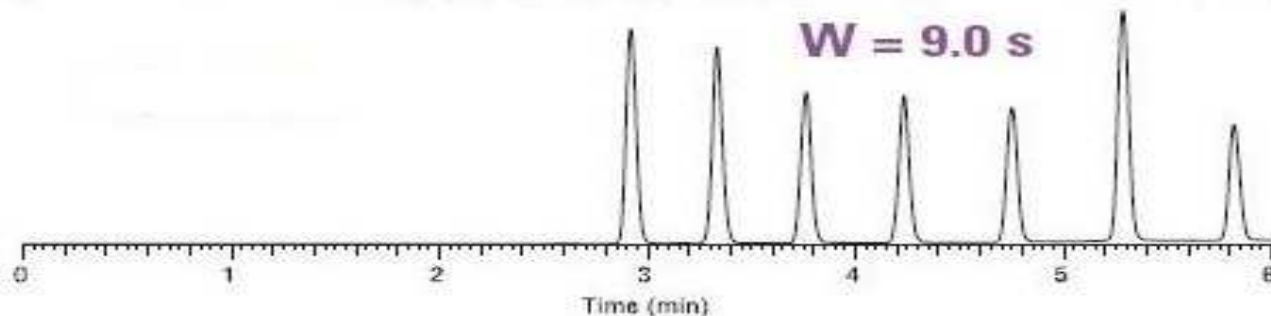
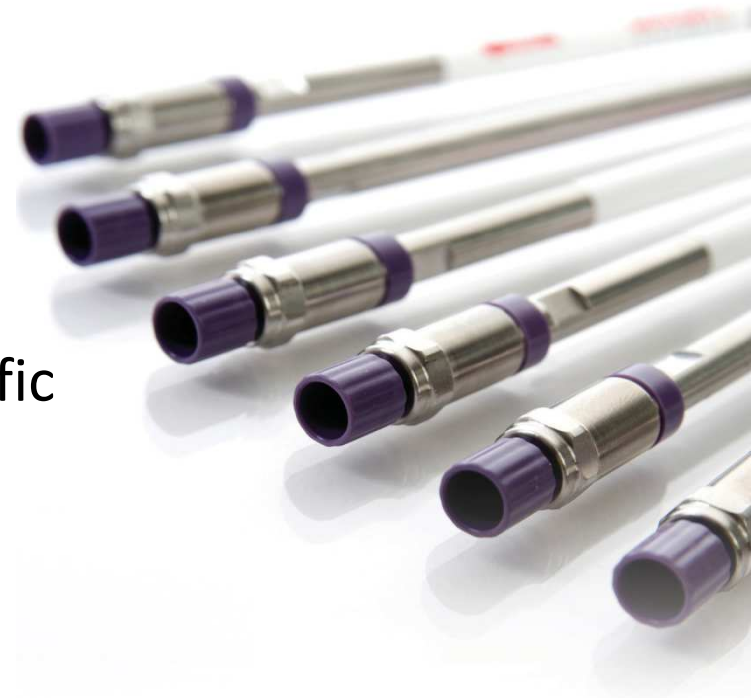
Comparisons

■ Existing method
 ■ New Column Dimensions
 ■ Adjust Flow Rate

Existing method vs. New Column Speed (unit) Pressure (unit)	Existing method vs. Variable Flow Speed (unit) Pressure (unit)	Time Saved Existing Method (minutes) New Column (minutes) Variable Flow (minutes)	Solvent Comparison Existing Method (ml) New Column (ml) Variable Flow (ml)
--	---	---	--

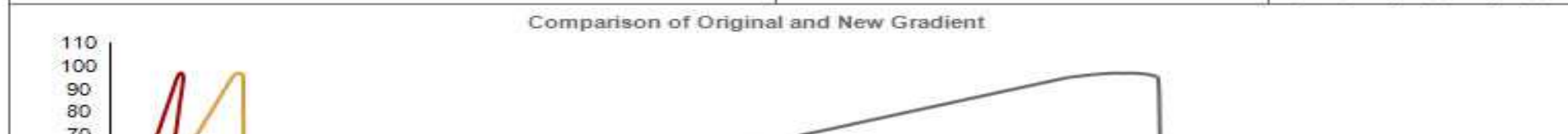Virtuoso Vial Identification System

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LC Columns and Accessories

Section Contents

Introduction	4-001	LC Accessories	4-158
Column Selection	4-002	LC Equipment	4-186
Column Protection	4-016	LC Reagents	4-193
Columns for Fast and Analytical LC	4-022	LC and LC-MS Instrument Parts	4-209
Columns for Biomolecules	4-110	Technical Information	4-231

Featured Products

Accucore Columns

Ultimate core performance - speed and selectivity combined

➤ [PAGE 4-042](#)

Synchronis Columns

Consistent, reproducible separations, column after column, time after time

➤ [PAGE 4-074](#)

Hypersil GOLD Columns

Outstanding peak shape for your separations

➤ [PAGE 4-024](#)

MABPac Columns

Unrivalled resolution and efficiency in the analysis of protein variants

➤ [PAGE 4-126](#)

Acclaim PepMap Columns

The standard for peptide separations in proteomics

➤ [PAGE 4-149](#)

Viper Connectors

Simple, dead volume free plumbing of HPLC and UHPLC systems

➤ [PAGE 4-158](#)

← **Table of Contents** ✕

- Introduction
- Column Selection
- Column Protection
- Columns for Fast and Analytical LC
- Application Specific LC Columns
- Columns for Biomolecules
- LC Accessories
- LC Equipment
- LC Reagents
- LC Instrument Parts
- Technical Information

Contents
Search

Section Contents

Introduction	4-001	LC Accessories	4-158
Column Selection	4-002	LC Equipment	4-186
Column Protection	4-016	LC Reagents	4-193
Columns for Fast and Analytical LC	4-022	LC and LC-MS Instrument Parts	4-209
Columns for Biomolecules	4-110	Technical Information	4-231

LC Columns and Accessories

Products

Synchronis Columns

Consistent, reproducible separations; column after column, time after time

➤➤ [PAGE 4-074](#)

Hypersil GOLD Columns

Outstanding peak shape for your separations

➤➤ [PAGE 4-024](#)

Thermo Scientific
Chromatography Columns
and Consumables 2014-2015

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