RHEOLASER CRYSTAL



 \bigcirc

MONITOR ANY PHYSICAL PHENOMENON Crystallisation, melting, polymorphic transition...

<u>^</u>

ACCURATE 4 – 90°C 0.1 – 25°C/min Optical non-invasive measurement

VERSATILE

 $\overline{}$

Any sample size (0.05-5g) Any sample structure Any sample form

$\widehat{}$

STRAIGHTFORWARD 1-click measurement Direct sampling No denaturation Disposable sample holders

MICROSTRUCTURE EVOLUTION THERMAL ANALYSIS BY MICRORHEOLOGY




MICRORHEOLOGY

www.formulaction.com

RHEOLASER CRYSTAL


TAKE A DEEPER LOOK AT YOUR SAMPLES MICROSTRUCTURE

Rheolaser Crystal enables to monitor microstructure evolution in heterogeneous products with an innovative optical method, combining a non-invasive measurement with accurate temperature control and sufficiently large sample volume (0.05 - 5g) to overcome problems of heterogeneities. This enables to measure finished products, such as food, cosmetics or pharmaceuticals, and identify transition temperatures of proteins, polymers, waxes or any fatty compounds.

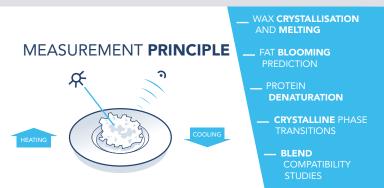

Rheolaser Crystal uses the Diffusing Wave Spectroscopy principle (DWS). Light is scattered by the particles, creating an interference pattern (Speckle Image). The variation of this image is directly related to the motion of the particles: the faster they move, the faster the Speckle Image changes. By a mathematical analysis of this variation, decorrelation functions can be computed and then processed, to obtain a characteristic time τ , as a function of time or temperature.

Values of $1/\tau$, or **Micro-Dynamics (Hz)**, are then plotted against time or temperature, resulting in characteristic peaks when the product shows a microstructural evolution, such as a phase transition or any other physical event. The signal can then be integrated for an easier visualisation, obtaining the so-called **Micro-Dynamics Evolution (%)**.

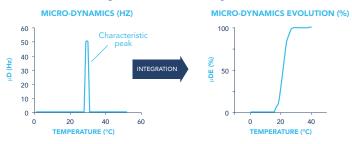

KEY **BENEFITS**

A VERSATILE DEVICE

Any sample can be analysed, no matter its form (liquid, solid, gel...) or the volume available (from 0.05g for precious products, to 5g for heterogeneous samples). Use the instrument for complex mixtures and blending compatibility studies.

CHARACTERISE YOUR SAMPLE AS IT IS

Thanks to the non-intrusive measurement, and to the large and direct sampling, there is absolutely no risk to damage the structure of the sample, or to obtain non-representative data because of heterogeneities. Rheolaser Crystal will just measure your whole sample.

FAST & ACCURATE ANALYSIS

Analysis can be launched in just 1-click, after direct sampling of the product. The optical DWS principle, enabling sensitivity at the nanometer size, combined to a very accurate temperature control, makes Rheolaser Crystal the fastest and most accurate way to monitor microstructure evolution.



TAL

APPLI**CATIONS**



TECHNICAL SPECIFICATIONS

Temperature range	4 – 90°C	
Temperature ramp speed	0.1 – 25°C/min	
Sample volume	0.05 – 5 g	
Sample form	Any (liquid, solid, gel)	
Disposable sample holder	•	
Laser wavelength	650 nm	
Dimensions (cm)	47x26x34	
Weight (kg)	15 kg	
Recommended configuration	Windows 7, 8 or 10, Intel Core i5, 4 Gb RAM, 32/64 bit	

