



HPLC columns

Acclaim HPLC columns

Product manual

Contents

Introduction	
Introduction to the Acclaim HPLC column	3
Systems and instrumentation requirements	3
Before you start	
The most important rules	4
Acclaim column operational parameters	4
Operation and system requirements	
General system requirements	5
Installation and start-up	
System rinse	6
Installing the column	6
The quality assurance report	6
Preparation of eluents and standards	
Mobile phase preparation	7
Sample preparation	7
Column care	8
Column storage-long term storage	8
Mobile phase	8
Recommended operating pH range	8
Recommended operating temperature limit	8
Flow rate and pressure limit	8
Cleaning the column	8
Ordering information	9

Introduction

Introduction to the Acclaim HPLC column

The Thermo Scientific™ Acclaim™ HPLC Columns are based on spherical, high-purity silica particles. The columns feature a well-balanced integration of high column efficiency, excellent performance, complementary selectivity, as well as stable and rugged column packing. The Acclaim columns packed with 3 µm silica are designed to generate less backpressure compared to other small particle columns, and are more resistant to column fouling, making them compatible with standard HPLC instrumentation. The Acclaim columns packed with 2.2 µm silica are designed for very high resolution separations and are compatible with ultra-high pressure instrumentation. As the result, Acclaim columns provide rapid separation solutions in a broad range of applications, including pharmaceutical, food and beverage, environmental, chemical, consumer products, and more.

The Acclaim columns are silica based, 2.2 µm and 3 µm particle size. There are four stationary phases: Acclaim 120 C18, Acclaim 120 C8, Acclaim Polar Advantage and Polar Advantage II. The Acclaim C18 and C8 columns (2.2 µm and 3 µm), feature high-density monomeric C18 and C8 chemistry (respectively) with exhaustive end-capping, and provide the same selectivity as the standard phases and extremely low silanol activity. Acclaim Polar Advantage column is a unique polar-embedded phase compatible with aqueous mobile phases containing no organic solvent modifiers, though maintaining selectivity similar to C18 column. Acclaim Polar Advantage II column features amide-embedded chemistry with excellent hydrolytic stability (pH 1.5 to 10), and provides complementary selectivity.

The manufacture of all Acclaim columns is controlled and documented according to ISO 9001 guidelines. Certificates of Quality Assurance and Lot Validation accompany each analytical column.

Systems and instrumentation requirements

The Acclaim HPLC columns allow accelerated separations to be run on standard LC instrumentation, provided these instruments fulfill distinct minimum requirements. In order to achieve the optimal speed of analysis and resolution, it is advantageous to use systems with lowest gradient delay volume and extra column volume. The detector must provide a minimum data rate and the flow cell needs to support these requirements for extra column volume. We suggest the use of one of Thermo Scientific™ HPLC/UHPLC systems that meets the requirements.



Before you start

The most important rules

This manual describes the recommended operating conditions for the Acclaim HPLC columns, and the precautions to maximize and preserve its performance.

1. Observe the recommended ranges of operating conditions.
2. Keep the column clean by using HPLC-grade filtered mobile phases, by filtering the samples, by using pre-column filters.
3. Pay close attention to minimizing dead volume in the various connections, and gradient delay volume in the pump
4. Equilibrate the column sufficiently before use, and store it properly after use.

Acclaim column operational parameters

Table 1. Recommended ranges of operation

Parameter	Acclaim 120 C18 column	Acclaim 120 C8 column	Acclaim Polar Advantage column	Acclaim Polar Advantage II column
pH range	2.0–8.0	2.0–8.0	2.0–8.0	1.5–10
Temperature	60 °C	60 °C	60 °C	60 °C
Aqueous compatibility	0–90%	0–90%	0–100%	0–100%

The type 316 stainless steel column hardware is compatible with common HPLC mobile phases. To prevent corrosion, high concentrations of halide ions at pH<2.5 should be avoided. The pressure limits are dependent on the column geometry and particle size, and are the same for all surface chemistries.

Table 2.

Particle size	Geometry	Recommended maximum pressure	Recommended maximum flow rate	Typical flow rate
3 µm	3.0 x 33 mm	4,500 psi	2.5 mL/min	0.4 – 1.0 mL/min
	3.0 x 50 mm	5,000 psi	2.5 mL/min	0.4 – 1.0 mL/min
	3.0 x 75 mm	6,000 psi	2.5 mL/min	0.4 – 1.0 mL/min
2.2 µm	3.0 x 30 mm	4,500 psi	2.5 mL/min	0.6 – 1.6 mL/min
	3.0 x 50 mm	7,000 psi	2.5 mL/min	0.6 – 1.6 mL/min
	3.0 x 100 mm	10,000 psi	2.5 mL/min	0.6 – 1.6 mL/min
	2.1 x 30 mm	4,500 psi	1.2 mL/min	0.3 – 0.8 mL/min
	2.1 x 50 mm	8,500 psi	1.2 mL/min	0.3 – 0.8 mL/min
	2.1 x 100 mm	11,000 psi	1.2 mL/min	0.3 – 0.8 mL/min
	2.1 x 150 mm	11,000 psi	1.2 mL/min	0.3 – 0.8 mL/min
	2.1 x 250 mm	15,000 psi	1.2 mL/min	0.3 – 0.8 mL/min

Operation and system requirements

General system requirements

The Acclaim HPLC columns can provide significant benefits to any standard HPLC system, but for the best results we recommend systems with low extra-column dispersion the following minimum requirements must be fulfilled:

- System gradient delay volume not larger than 1200 μL .
- Connection tubing internal diameter not wider than 250 μm (0.010 inch).
- Flow rates up to 3 mL/min for long term use.
- Upper pressure limit of at least 350 bar (5000 psi).
- Column thermostating ability with minimum extra column contribution.
- Detector flow cell volume not larger than 13 μL .
- Detector data collection rate of minimum 10 Hz.
- Detector time constant must be settable to values of as low as 0.2 seconds.

Pump

For isocratic methods, the choice of pump is not critical. For gradient elution, the gradient delay volume is a significant factor, and should be minimized. Gradient delay volume is the volume between where the solvents first mix to the head of the column. The low-pressure gradient-mixing (LPG) design is economical and flexible, but has a greater delay volume; the high-pressure gradient-mixing (HPG) design while more expensive has much lower gradient delay volume. Gradient mixing devices can make a large contribution to the delay volume, but are usually necessary for best results; therefore choose the smallest one that is consistent with good gradient mixing in the flow range of 0.4 to 1.3 mL/min.

Injector

When moving to a 3.0 mm ID column from 4.6 mm, you normally reduce the injection volume by a factor of 0.4. The popular split-loop autosampler design puts a relatively large sample loop in line; to reduce gradient delay volume, you may want to install a smaller loop, or program the valve to bypass the loop shortly after injection.

Detector

The flow cell makes little difference in peak shape when the peak volume is greater than 10 $\times$ the flow cell volume. Since the peak volume (width at base times flow rate) is reduced to as little as 20% of the original when you accelerate a method using 2.2 μm particles, you should consider if a smaller flow cell is appropriate. For good integration of peaks, you should have 20 – 30 data points across the narrowest peak of interest, therefore set the data collection rate and filtering constant appropriately.

Tubing connections

To minimize extra-column volume, use the recommended size or smaller capillary connections between the column and injector, and between the column and detector. Make sure that the tubing is cut squarely and cleanly, and that all fittings are correctly made. The use of Thermo Scientific™ Viper™ Fingertight Fitting Systems is strongly recommended.

Column thermostat

This module is optional but recommended. If you are operating more than 15 °C different than ambient, you should consider a mobile phase pre-heater.

Pre-column filter

Guard columns are designed to protect the main analytical column in chromatography by trapping contaminants and particulates from the sample. This is especially useful when working with dirty or complex samples.

Installation and start-up

System rinse

Prepare fresh, clean mobile phases. Prime the pump with the mobile phases to make sure that all air and previous mobile phases are expelled.

Installing the column

1. The columns are shipped in an acetonitrile-water mixture, therefore be sure that the mobile phase is compatible with that. Before installing the column, purge all the pump lines with liquid to remove air and ensure the system is clean.
2. Connect the column between the injection valve and the detector, with the flow arrow on the label pointing towards the detector. The column end fittings are compatible with standard 10-32 fittings and ferrule dimensions. Use the shortest lengths of tubing practical to minimize system volume. The tubing ID should not exceed 0.007 inches (0.18 mm) for 3.0 mm ID columns nor exceed 0.005 inches (0.125 mm) for 2.1 mm columns. Thermo Scientific Viper Fingertight Fitting Systems are strongly recommended.
3. Once the column is installed, pump at least 10 column volumes of mobile phase through the column, and if additives such as ion pairing reagents are present in the mobile phase, up to 200 column volumes may be needed. The column is equilibrated when the baseline is stable and several injections produce stable retention times.

The quality assurance report

Test the Acclaim column prior to its first use. Using the conditions described on the Quality Assurance Certificate, confirm that it performs as indicated. Keep a record of its performance for future reference.

Preparation of eluents and standards

Mobile phase preparation

High purity chemicals

For best results use high-purity deionized water and HPLC-grade organic solvents; these will have the least UV background, and are pre-filtered. The Acclaim columns may be used with all the usual HPLC mobile phases.

Filtration

Buffer salts used for the preparation of mobile phases often contain insoluble particulates. If these are allowed to flow into the column, they will be trapped at the head of the column by the inlet frit, plugging the column over time. To avoid this the mobile phase should be filtered through a 0.5 µm or finer porosity filter media. In order to avoid problems associated with the loss of prime in the pump, mobile phases should be degassed prior to use.

Solvent compatibility

The columns are compatible with all commonly used HPLC solvents, including methanol, acetonitrile, tetrahydrofuran, acetone and dichloromethane. Acclaim 120 C18 and C8 columns are compatible with highly aqueous mobile phases (up to 90% aqueous) under isocratic conditions. In gradient methods, they are compatible with up to 95% aqueous mobile phases.

Thermo Fisher Scientific recommends Acclaim Polar Advantage and Polar AdvantageII columns for applications that require highly aqueous mobile phases (> 90% aqueous). For control of mobile phase pH between pH 2 and pH 7, dilute acids and buffer salts such as trifluoroacetate, phosphate, formate, acetate or TRIS may be used. When making alkaline buffers for Acclaim Polar Advantage II column, Thermo Fisher Scientific recommends amine-type buffers such as TRIS, ammonia or diethylamine instead of phosphate buffers. When adding organic solvent to salt and buffer solutions, it is important to remember this may result in a precipitation of the salt, often in the column, leading to possible damage to the column. Always check the solubility of buffer salts in solvent-buffer mixtures before running new methods. For very fast gradient elution, greater than 40% per minute, we recommend acetonitrile as the organic solvent because of its lower viscosity.

Sample preparation

Contamination of the column by particulate matter in the sample is a leading cause of column failure. Symptoms may include high pressure, poor symmetry, ghost peaks, or low efficiency. Pass the sample through a 0.5 µm or smaller porosity filter before injection. Biological samples are especially troublesome due to dissolved proteins that may precipitate some time after the initial preparation. Short columns are more sensitive to overload conditions than long columns. If the sample is dissolved in a solvent stronger than the mobile phase the chromatogram will be susceptible to "volume overload." If possible, reduce the solvent content in the prepared sample to no more than the initial composition of the mobile phase.

Column care

Column storage-long term storage

If the column will not be used for a week or more, flush the column with 10 column volumes of unbuffered mobile phase containing at least 70% organic solvent. Do not store it with acids, salts or buffers inside. Plug the column to keep the bed moist. Do not allow the column to dry out.

Mobile phase

Mobile phases should be freshly prepared every time. All chemicals and solvents should be at the highest available quality. All mobile phases should be filtered before use. In-line filters are recommended.

Recommended operating pH range

To obtain better column lifetime, it is highly recommended to use "silica friendly" mobile phases. The pH limits of the columns are listed in Table 1.

Recommended operating temperature limit

The recommended temperature limits of the columns are listed in Table 1, and depend on the nature of the bonded phase. Higher temperatures accelerate the loss of bonded phase and dissolution of the silica substrate.

Flow rate and pressure limit

To protect the column, observe the recommended maximum flow and pressure. Best chromatographic results are usually at more moderate flows. See Table 2 for details. It is extremely important not to impose sudden column pressure surge. Therefore increase flow rate gradually up to the desired flow rate.

Cleaning the column

Hydrophobic compounds

If the column becomes contaminated with hydrophobic compounds, it may be safely washed with strong organic solvents. Flush out any buffer salts with unbuffered mobile phase. For a more detailed washing procedure, please refer to the [column care guide](#).

Ion-pairing agents usually can be removed by washing with 80:20 methanol:water acidified with 0.1% formic acid. Quaternary amine agents may require extensive washing to restore the original column behavior, and it is preferable to dedicate the column to this type of mobile phase before proceeding with the recommended wash procedure.

Particulate contaminants will cause excess backpressure and possibly loss of performance. Temporarily reverse the column, and at a reduced flow rate direct the effluent to waste until the pressure stabilizes. Prevention is the best cure: filter the samples and mobile phases, and use a pre-column filter.

Metal contamination (most commonly iron) may be removed by treatment with a chelating agent. The least damaging agent is 85:10:5 methanol:water:pentanedione. A more aggressive agent is 10 mM citrate, 20 mM EDTA at pH 4 mixed 70:30 with methanol.

Note: if above treatment fails to improve the column performance, replace it with a new one.

Ordering information

Acclaim HPLC columns ordering information

Column chemistry	Particle size	Column format	Quantity	Cat. no
120 C18	2.2 µm	2.1 x 30 mm	Each	071400
		2.1 x 50 mm	Each	068981
		2.1 x 100 mm	Each	068982
		2.1 x 150 mm	Each	071399
		2.1 x 250 mm	Each	074812
		3.0 x 50 mm	Each	071605
		3.0 x 100 mm	Each	071604
	3 µm	3.0 x 33 mm	Each	066272
		3.0 x 75 mm	Each	066273
120 C8	2.2 µm	2.1 x 100 mm	Each	072616
		2.1 x 150 mm	Each	072617
Polar Advantage	2.2 µm	2.1 x 50 mm	Each	072622
		2.1 x 100 mm	Each	072623
		2.1 x 150 mm	Each	072624
Polar Advantage II	2.2 µm	2.1 x 30 mm	Each	071402
		2.1 x 50 mm	Each	068989
		2.1 x 100 mm	Each	068990
		2.1 x 150 mm	Each	071401
		2.1 x 250 mm	Each	074814
		3.0 x 100 mm	Each	071607
	3 µm	3.0 x 33 mm	Each	066276
		3.0 x 50 mm	Each	068973
		3.0 x 75 mm	Each	066277

Guards and holder ordering information

For	Particle size	Column format	Quantity	Cat. no
Acclaim 120 C18 column	5 µm	3.0 x 10 mm	2/pack	071981
Acclaim 120 C8 column	5 µm	3.0 x 10 mm	2/pack	071979
Acclaim Polar Advantage column	5 µm	3.0 x 10 mm	2/pack	071983
Acclaim Polar Advantage II column	5 µm	3.0 x 10 mm	2/pack	071985
Acclaim Holder V-2 for 3 mm guard	–	–	Each	069580

 Learn more at thermofisher.com/acclaim