

7th Czech Mass Spectrometry
Conference

April 11th, 2018

ThermoFisher
SCIENTIFIC


Next Generation Technology for Reproducible and Precise Proteome Profiling

Lars Kristensen, Ph.D.
Application and training specialist

The world leader in serving science


- 1) Introduction and Q Exactive HF-X**
- 2) High Resolution DIA Workflow**
- 3) DDA+ Workflow (Label free quantification)**

Large Scale Proteomics

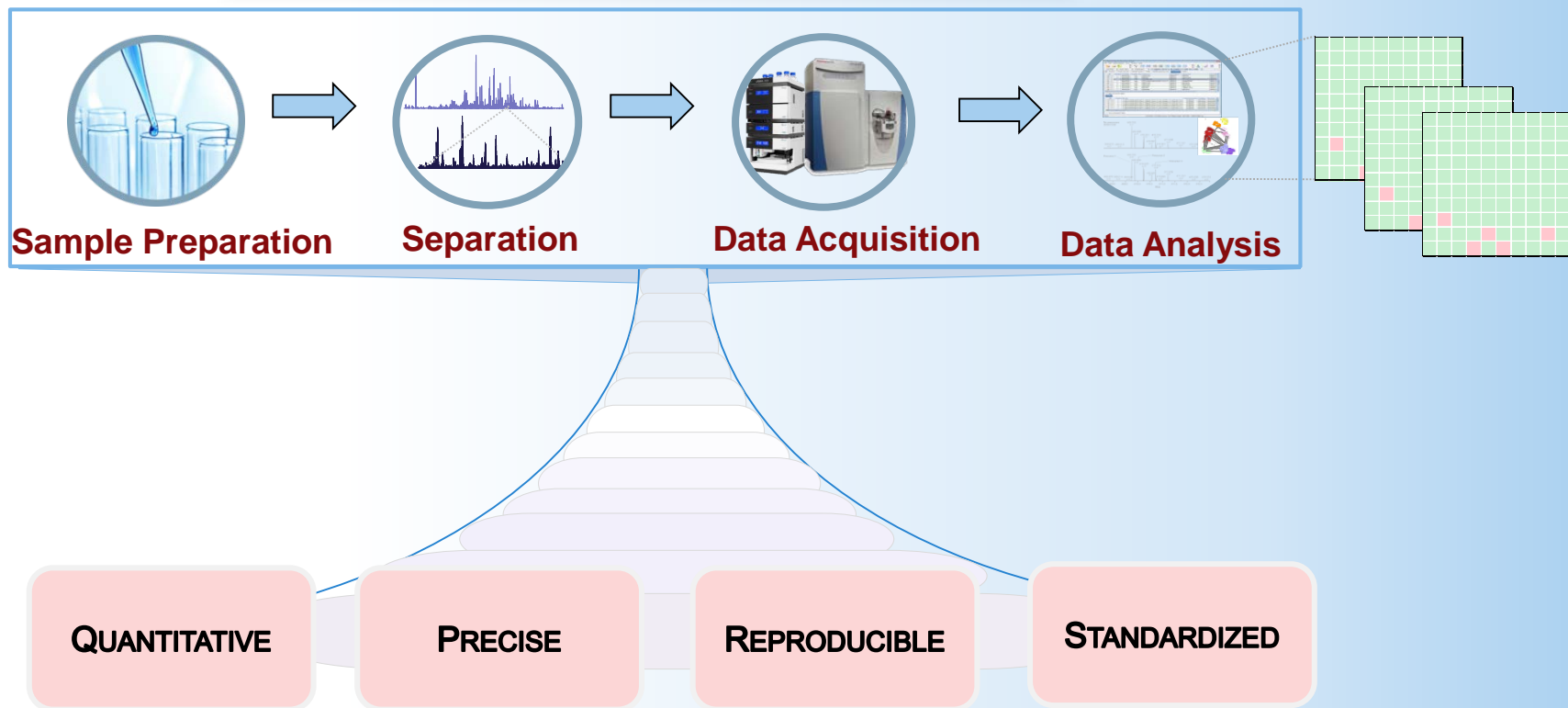

High-Resolution DIA Workflow

Unparalleled proteome coverage and dynamic range

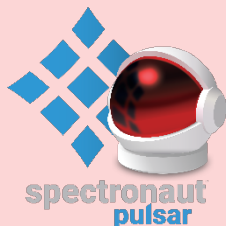

- Highest depth of proteome coverage and quantitative insight
- Robust quantitative precision

- ***Biospecimen profiling***
- ***Digital archiving***

DDA+ Workflow

Unsurpassed quantitative precision and reproducibility



- Unrivaled precision in precursor quantitation
- Maximize complete, reproducible quantitation across samples
- Minimize 'missing values' among samples

- ***Cellular signaling studies***
- ***Mechanism of action studies***
- ***PTM profiling***

Q Exactive HF-X – New architecture

Optimized Scan Matrix with accelerated HCD: 40 Hz MS/MS

Advanced DDA for bottom-up and top-down: Advanced Peak Determination

HyperQuad Mass Filter with Advanced Quadrupole Technology

Advanced Active Beam Guide

Electrodynamic Ion Funnel

C-Trap

HCD Cell


Ultra-High Field Orbitrap Mass Analyzer

High Capacity Transfer Tube

Dedicated transient for TMT 10plex: 50 K res


Improved S/N for intact proteins: 7.5 K res

Brighter ion beam: 2-3x More sensitive

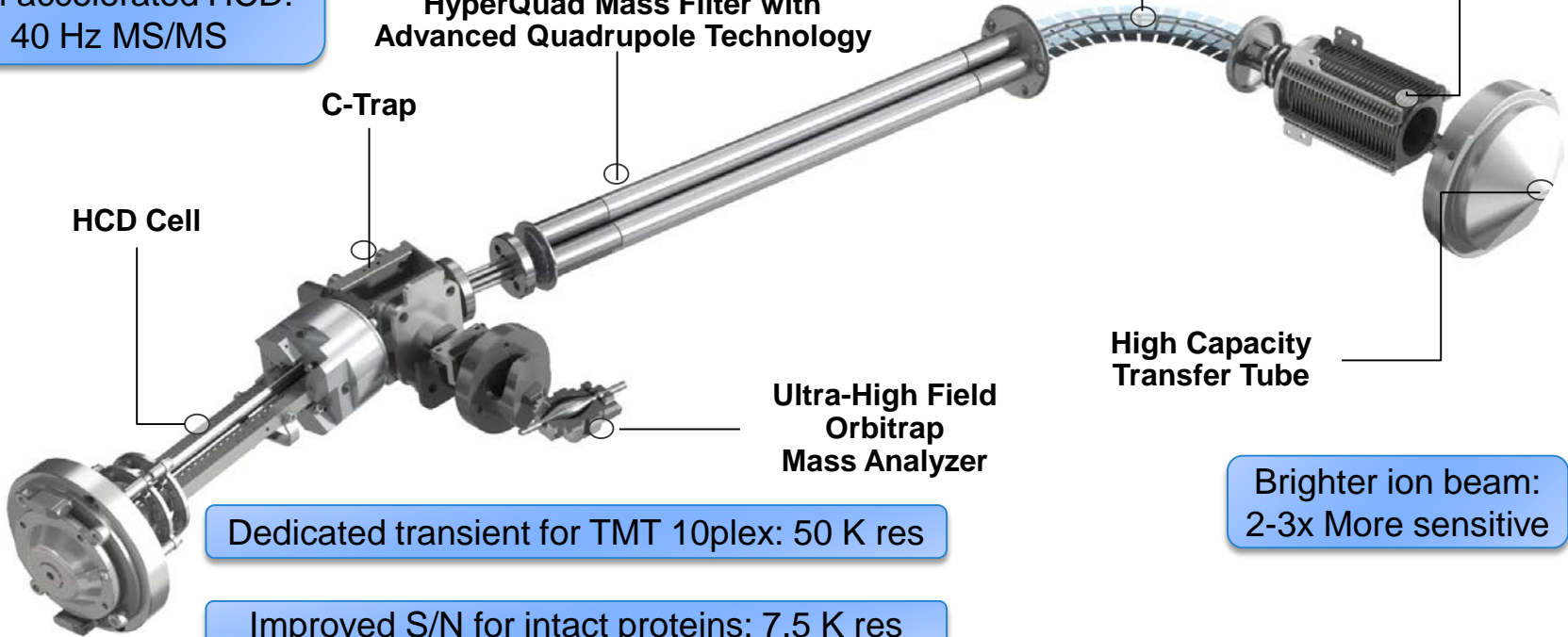

Q Exactive HF-X – Basis for increased speed and sensitivity


Q Exactive HF

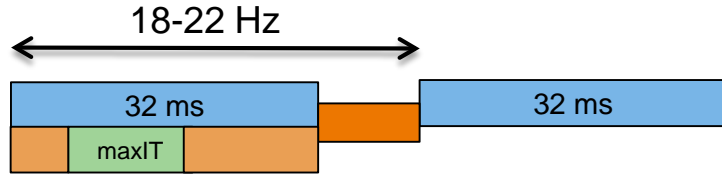


MS/MS

Longer fill time
Reduced scan overhead
aHCD


Q Exactive HF-X

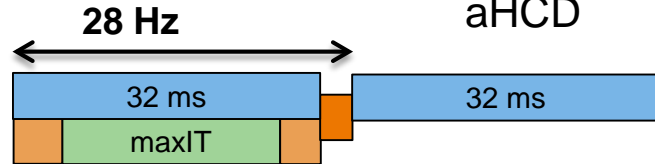


Comparable fill time
Reduced scan overhead
aHDC

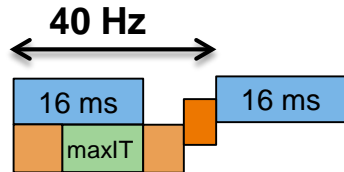


- Brighter ion beam (ion funnel), reduced scan overhead, and accelerated HCD (aHCD) is boosting acquisition speed
- Advantage for both MS and MS/MS mode
- Fast and high quality MS/MS acquisition up to 40 Hz with new 16 msec transient (7,500 resolution setting)

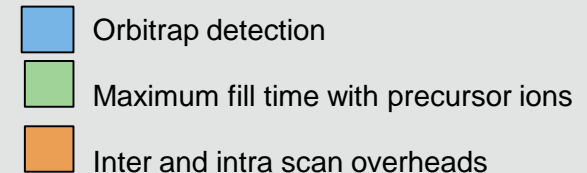

High-Resolution DIA Workflow

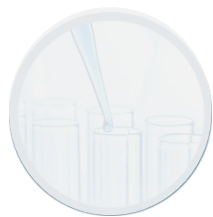


Sample Preparation

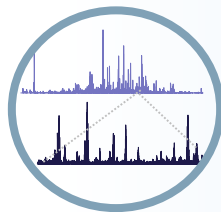


Separation



Data Acquisition

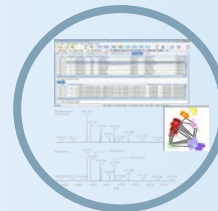

Data Analysis

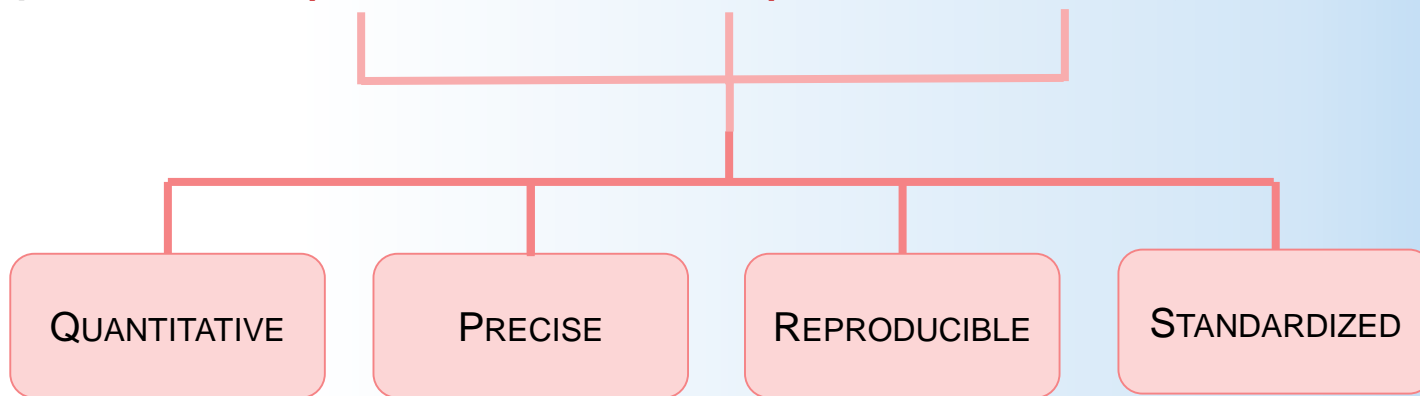

QUANTITATIVE

PRECISE

REPRODUCIBLE

STANDARDIZED

Workflow




Thermo Scientific™
UHPLC Systems



Thermo Scientific™
EASY-Spray™ LC Column



Thermo Scientific™ Q Exactive™ HF-X
Hybrid Quadrupole-Orbitrap MS

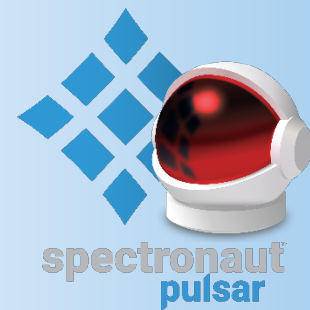


Spectronaut Pulsar™
software

- *Thermo Scientific™ UltiMate™ 3000 RSLCnano system*
- *Direct inject or pre-concentration mode*
- *Thermo Scientific™ Viper™ fittings*
- *150 μm ID x 150 mm,*
- *Sensitivity and robustness using Cap flow ($\approx 1.2 \mu\text{L}/\text{min}$)*
- *RT stability <1% observed for 350 injections*
- *Increased acquisition speed*
- *Advanced precursor determination*
- *Same # of protein IDs half the time*

Designed for Speed and Coverage

Key Benefits

- Spectronaut Pulsar™ is **specifically developed for the analysis of DIA & SWATH** data sets
- Data analysis with retention time correction based on spiked reference peptides using iRT Kit
- Spectral library generation from MaxQuant, Protein Pilot, and Thermo Scientific™ Proteome Discoverer™
- Direct visualization of qualitative and quantitative results on protein level including FDR and P-values
- Processing of very **large experiments** (1000s of runs)
- Fast data analysis speed in **less than 2 min per run**
- **DirectDIA™** library generation from DIA data (no need to create a DDA-based library)

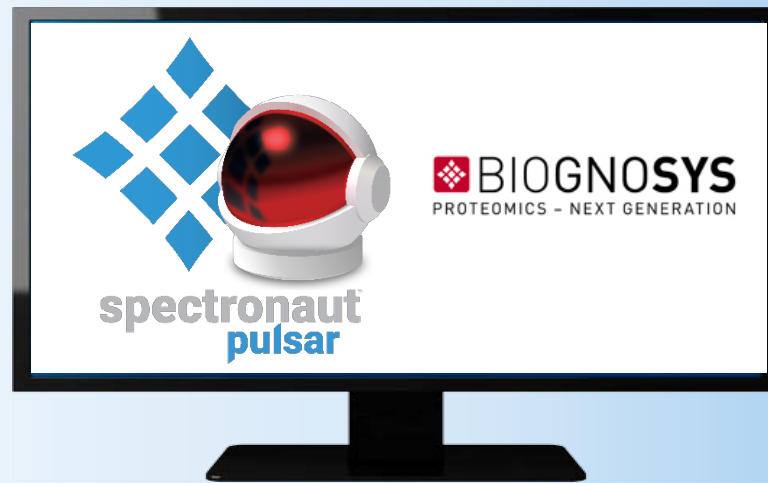

Designed for high throughput DIA data analysis





Balancing Efficiency Without Sacrificing Performance

Nanoflow

- Greater # of proteins
- Greater # of peptides
- Greater sensitivity
- Longer total run times


CapLC

- Greater Efficiency
- Shorter total run time (2X)
- Greater throughput
- More robust
- Less protein and peptide id's

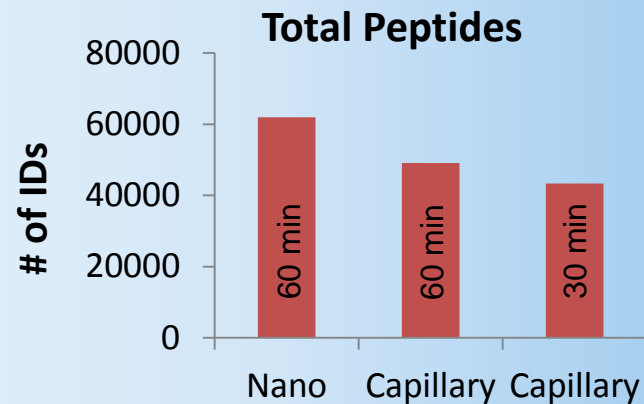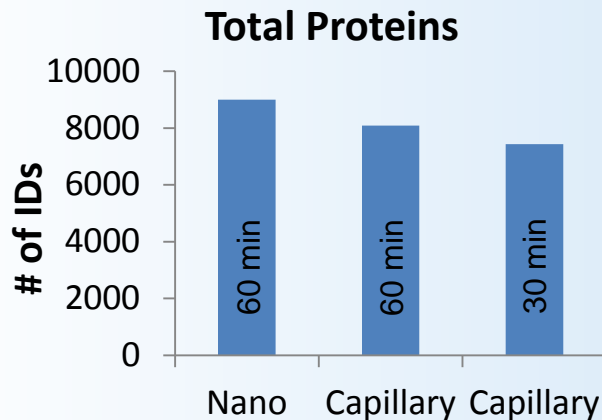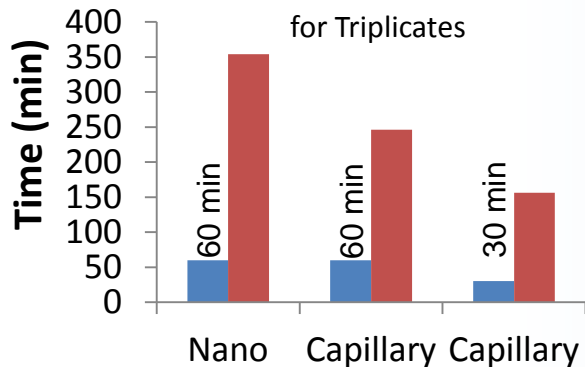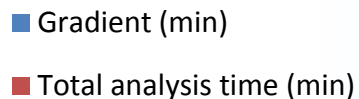


High-Resolution DIA Workflow: Highly Precise Proteome Quantitation


- Maximize depth of coverage
- Robust quantitative precision
- Confident in IDs
- Short analysis time

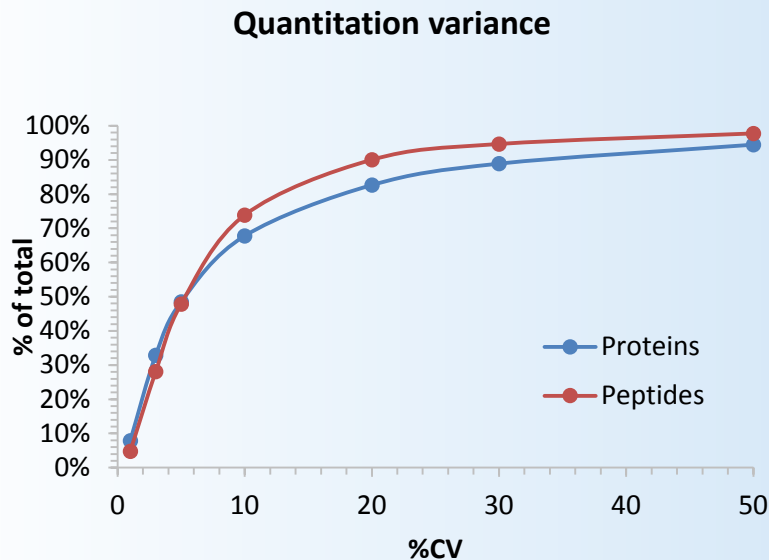


CapLC DIA, 4ug HeLa, 60min, 120K -> Spectronaut Analysis

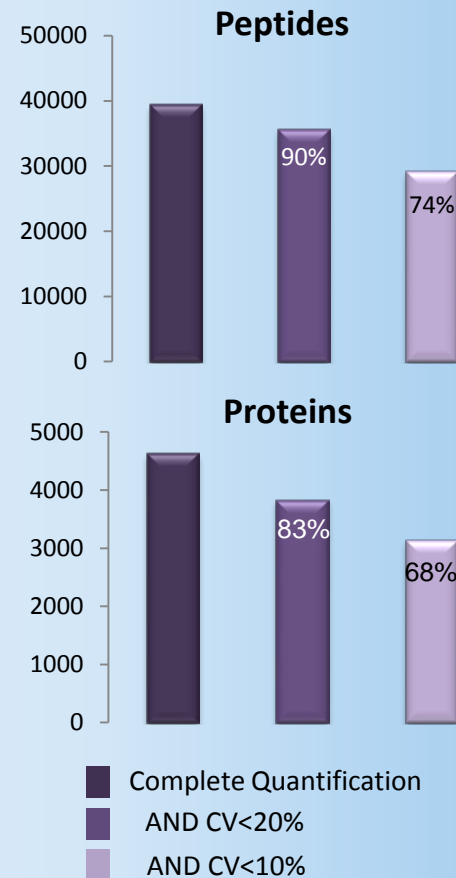

High-Resolution DDA+ Workflow

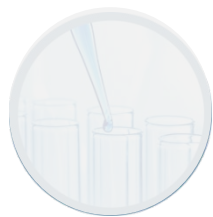


Sample Preparation

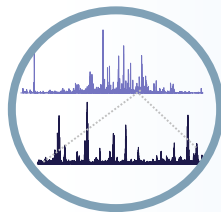


Separation



Data Acquisition

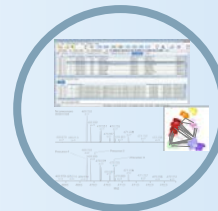

Data Analysis

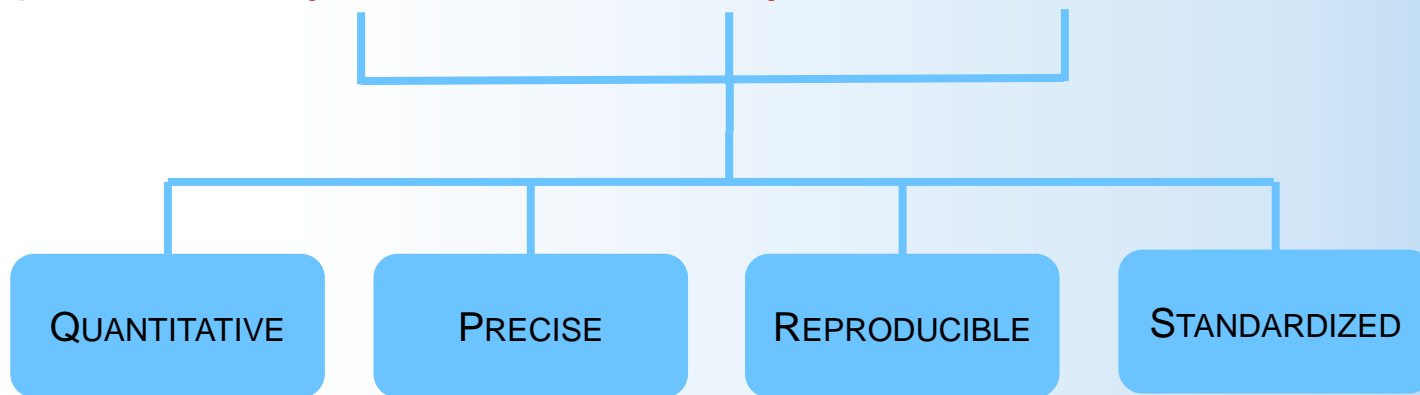

Workflow



UHPLC Systems

- *UltiMate 3000 RSLCnano system*
- *Direct inject or pre-concentration mode*
- *Viper fittings*



EASY-Spray LC Column

- *150 μm ID x 150 mm,*
- *Sensitivity and robustness using Cap flow ($\approx 1.2 \mu\text{L}/\text{min}$)*
- *RT stability <1% observed for 350 injections*



Q Exactive HF-X MS

- *Increased acquisition speed*
- *Advanced precursor determination*
- *Same # of protein IDs half the time*




Thermo Scientific™ Proteome Discoverer™ 2.2 software


Designed for Precision and Reproducibility

Key Benefits

- Enables large scale, multiplex proteomic studies (TMT 11-plex) and captures confident protein results which enables confident reproducibility
- Improved Label-free Quantitation
 - Feature mapping
 - Retention time alignment
 - Feature linking across files
- Minora Feature Detector node
 - Detects chromatographic peaks and features according to the specified quantification approach
- Minimizes 'missing data points' and maximizes quantitative insights

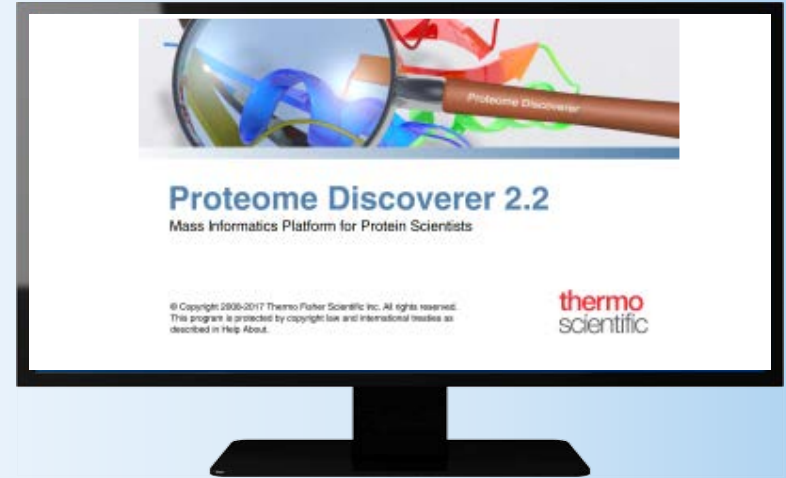


**Most comprehensive data analysis
platform for qualitative and
quantitative proteomics research**

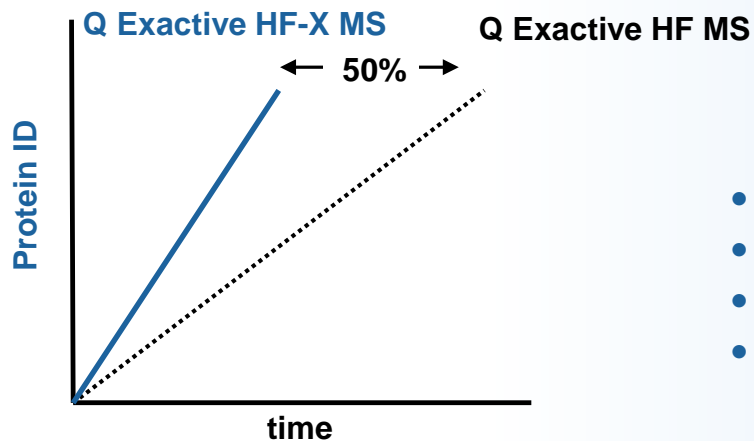

Maximizing protein identifications

- Quick screening of complex samples
- Quality control of complex samples
- Assessment of sample concentration
- Same # of protein ID in half the time

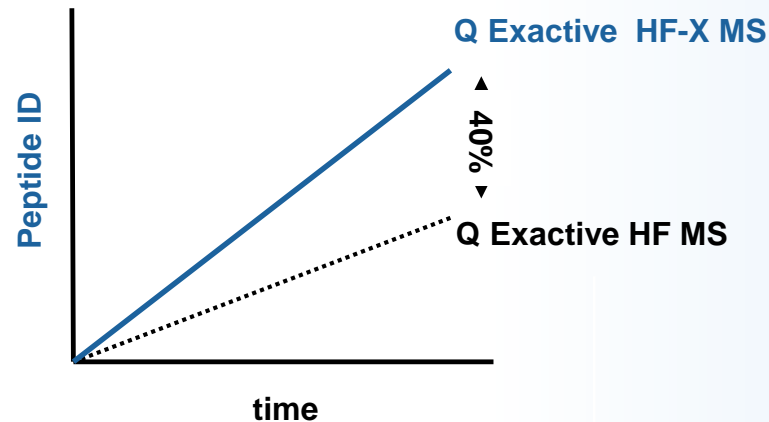


Maximizing peptide identifications

- Highest peptide coverage
- Deep proteome analysis
- Spectral library building

Saves time and samples in large-scale proteomics efforts

new

Completely quantified proteins

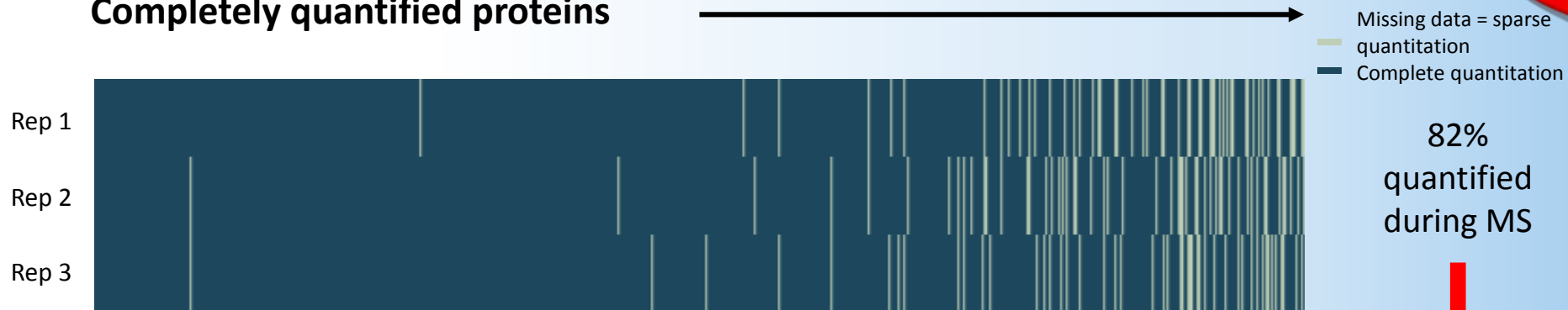

82% quantified during MS


↓

97% quantified from MS and Protein Discoverer 2.2 software

capLC DDA+, 4ug HeLa, 60min, 120K/7.5K, 19ms, Top 40 -> PD 2.2 Label-free Quant



Completely quantified peptides

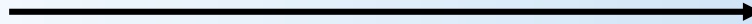


Missing data = sparse
— quantitation
— Complete quantitation

Rep 1

Rep 2

Rep 3

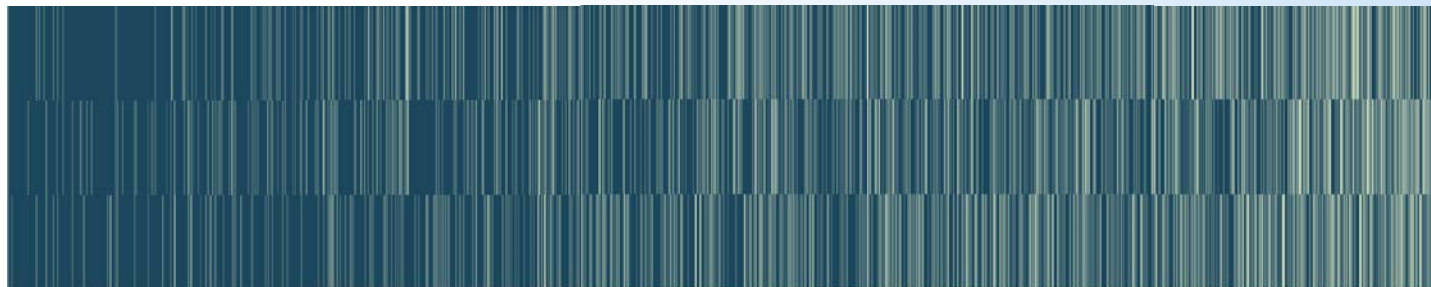

50%
quantified
during MS



97%
quantified
from MS and
PD 2.2

Rep 1

Rep 2

Rep 3



capLC DDA+, 4ug HeLa, 60min, 120K/7.5K, 19ms, Top 40 -> PD 2.2 Label-free Quant

Quantitation


Precision

Reproducibility


Standardization

DDA+ Enable Unrivalled Quantitative Precision

Protein quantitation variance

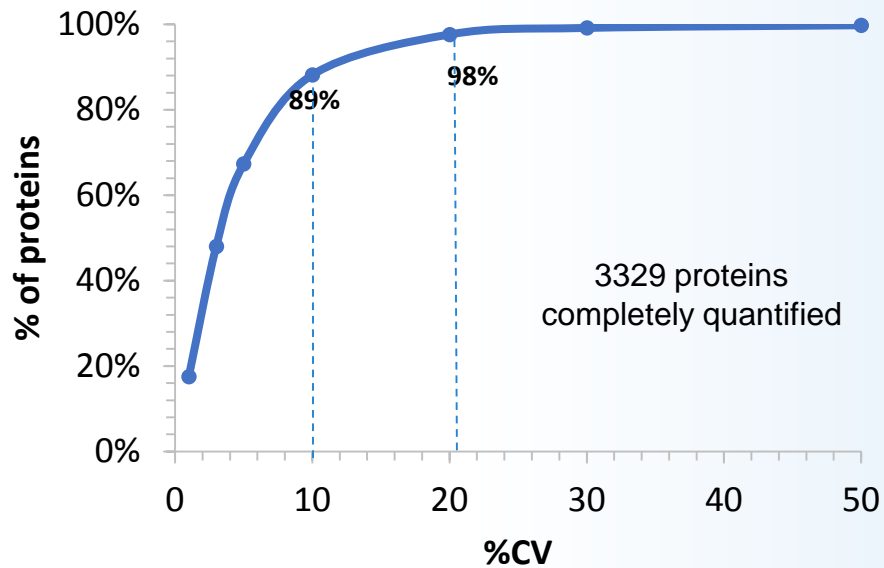

Peptide quantitation variance

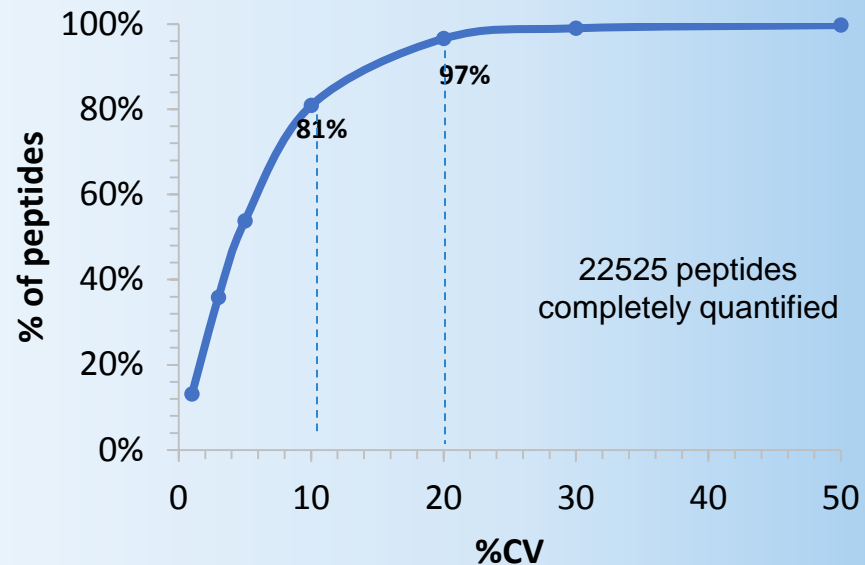

Quantitation


Precision

Reproducibility


Standardization

DDA+ Workflow: Greater Reproducibility Between Samples

Proteins

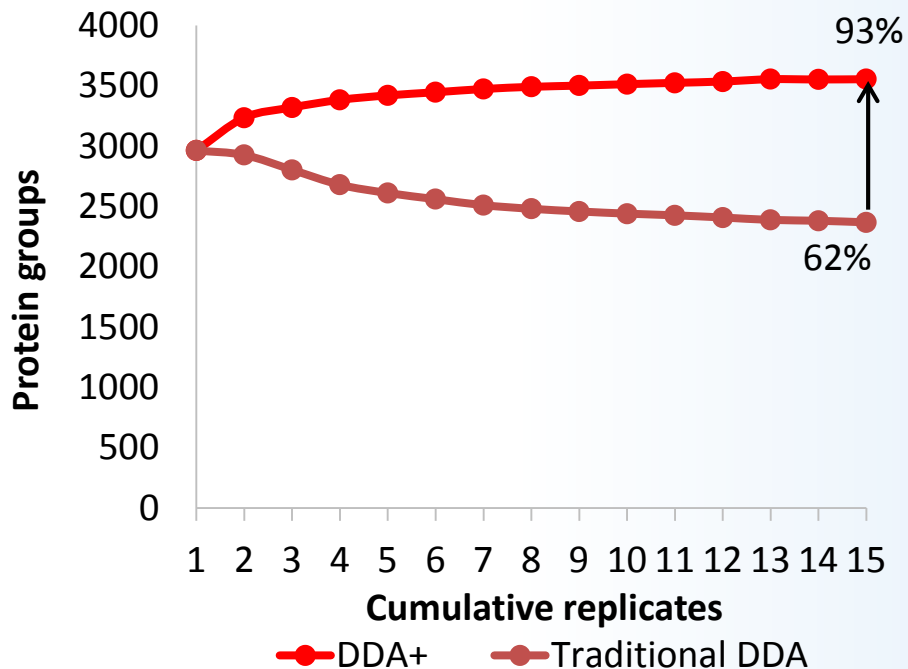

Peptides

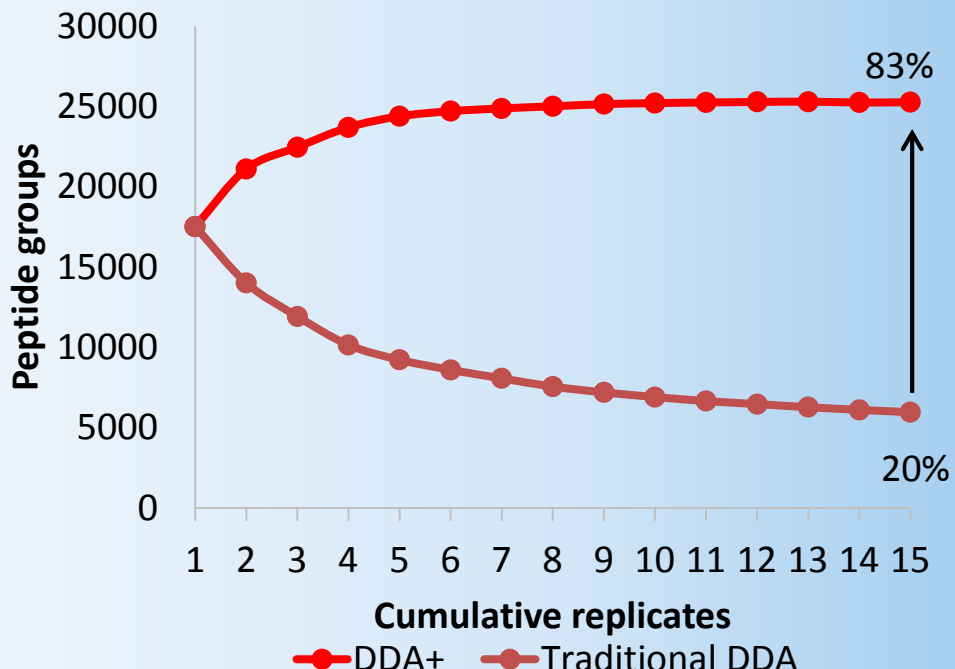

Quantitation

Precision


Reproducibility

Standardization

Inter-Site Consistency Across Different Instruments

Instrument Standardization Test

- Three Locations
- HeLa digest metrics
 - Protein
 - Peptide
 - PSMs
 - MS/MS
- 60 min gradient

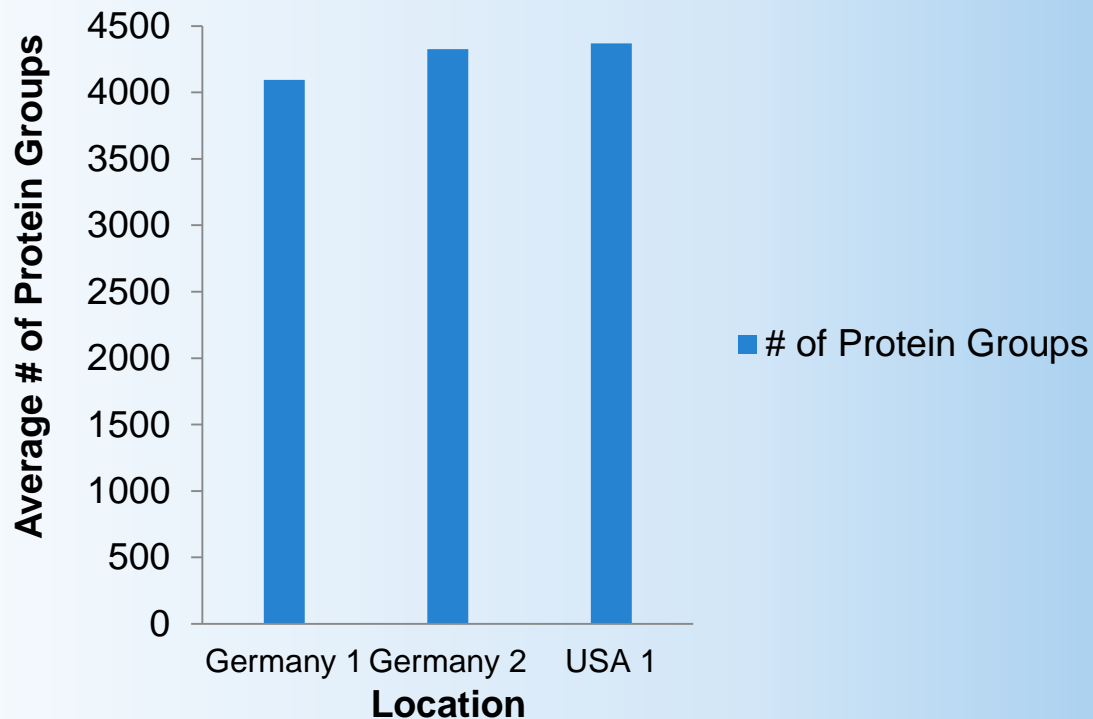



Questions?