pragolab

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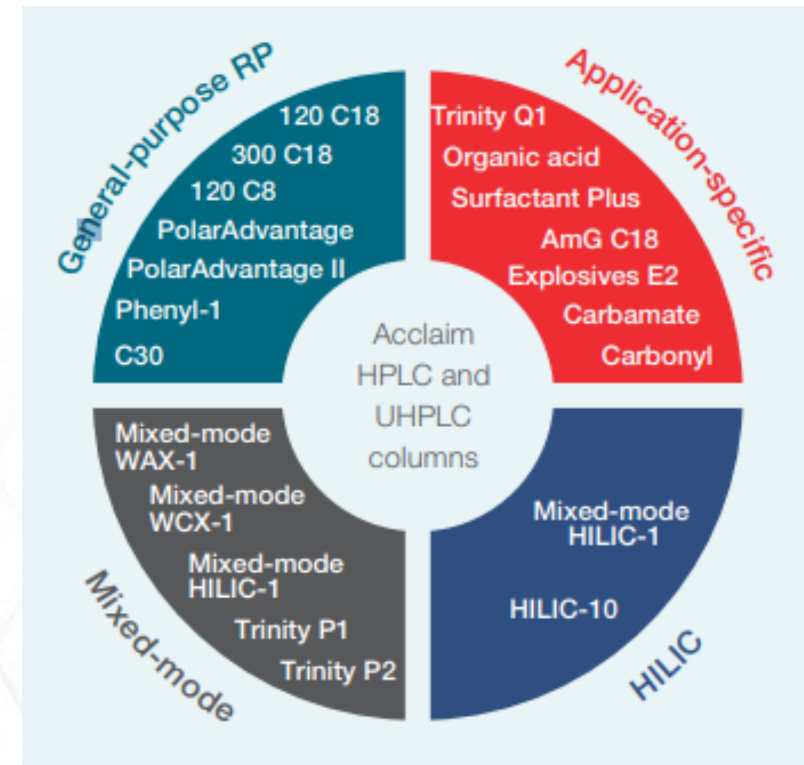
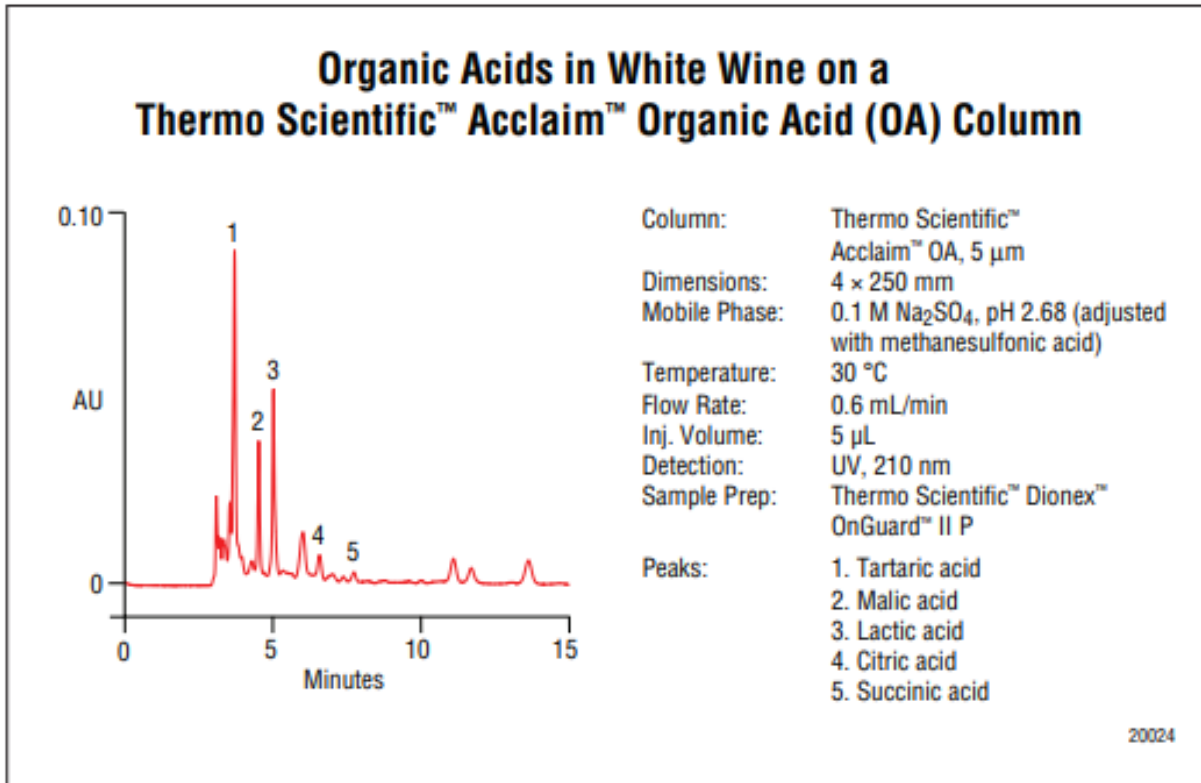
Contact Us

Sign In ▾

Welcome to the Thermo Scientific
AppsLab Library of Analytical Applications!

Find methods for your needs

enter compound, matrix or instrument type

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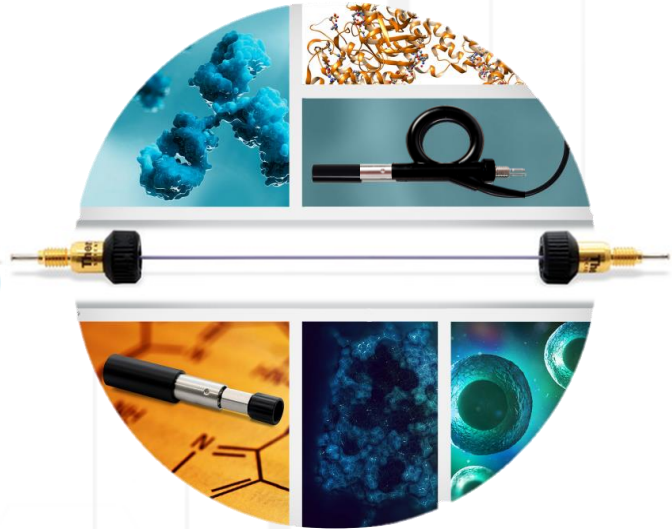
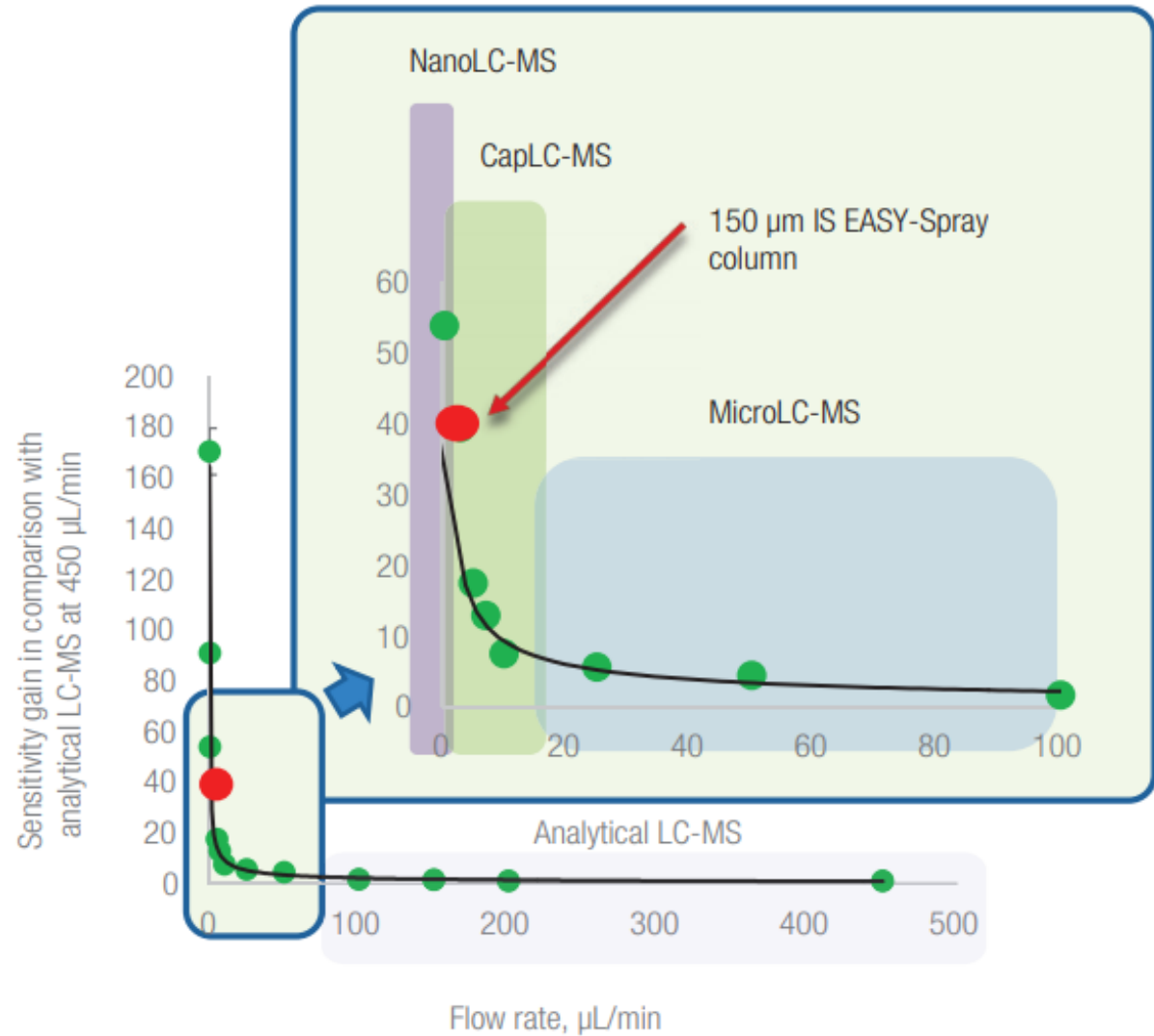
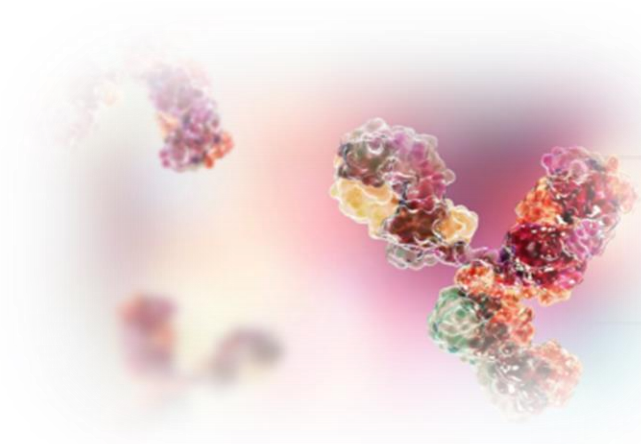
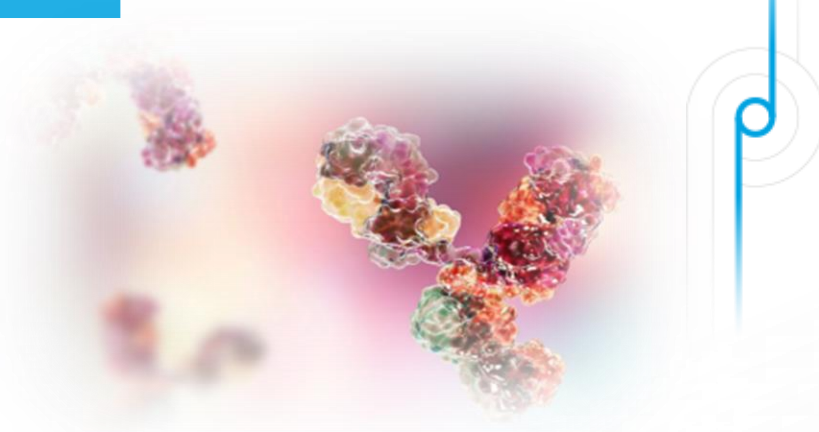
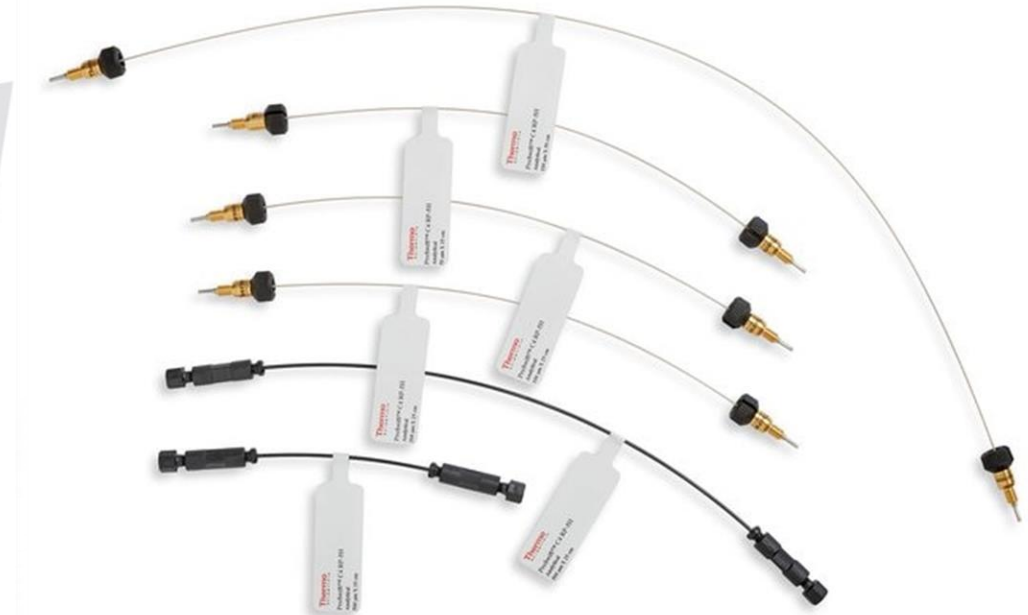
BIOCOMPATIBILITY

Available in biocompatible materials

MORE OPTIONS

A variety of materials and dimensions (length and ID) available for nearly every system and application

HIGH TEMPERATURE

Withstands temperatures up to 120 °C

REMOVABLE KNURL

Easy to grip and easy to remove, taking up less space to fit the tightest gaps

ID COLOR CODE

Easy visualization with color-coded tags for quick ID recognition

FINGERTIGHT

Easy installation and fingertight up to 1500 bar

IDENTIFICATION TAG

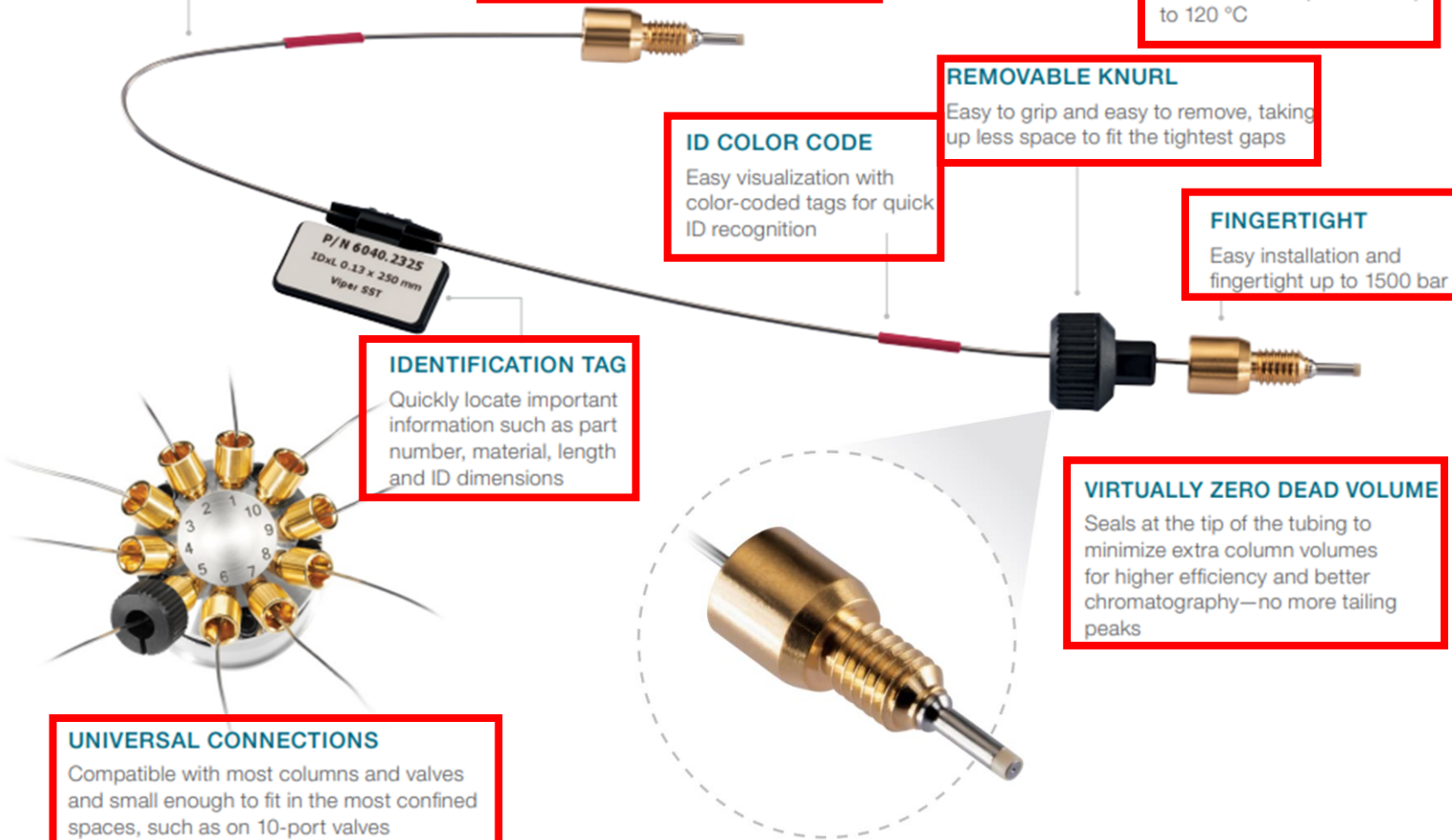
Quickly locate important information such as part number, material, length and ID dimensions

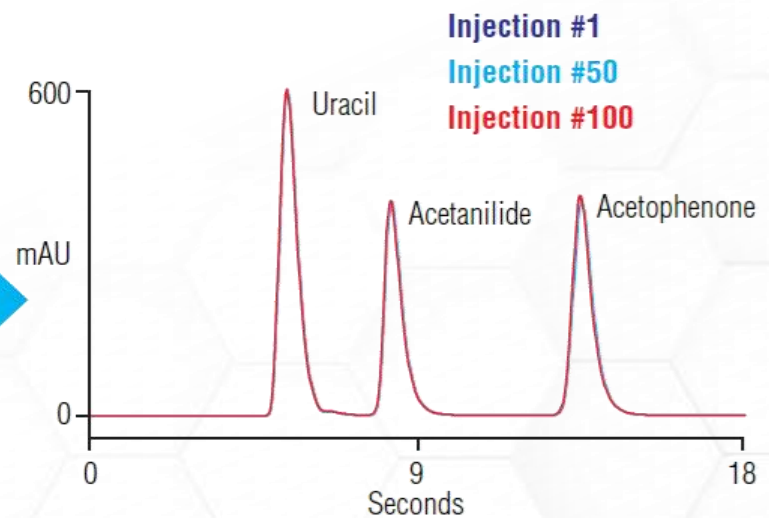
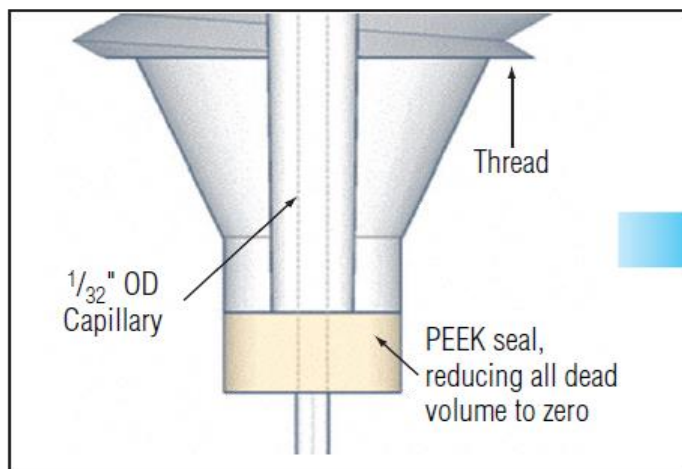
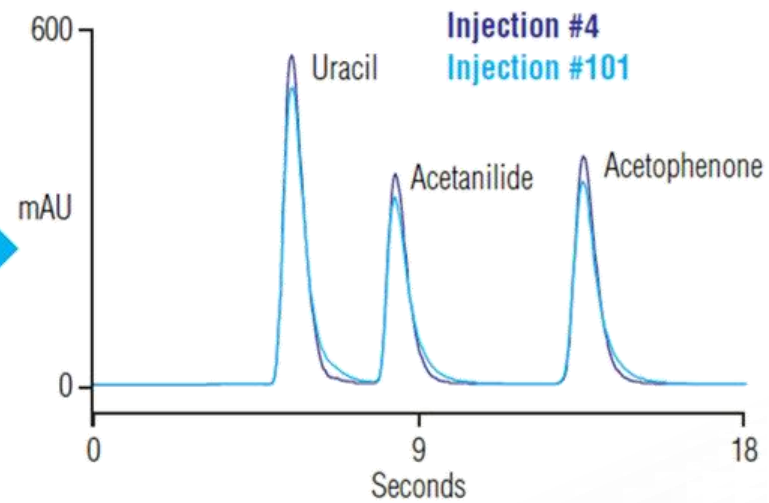
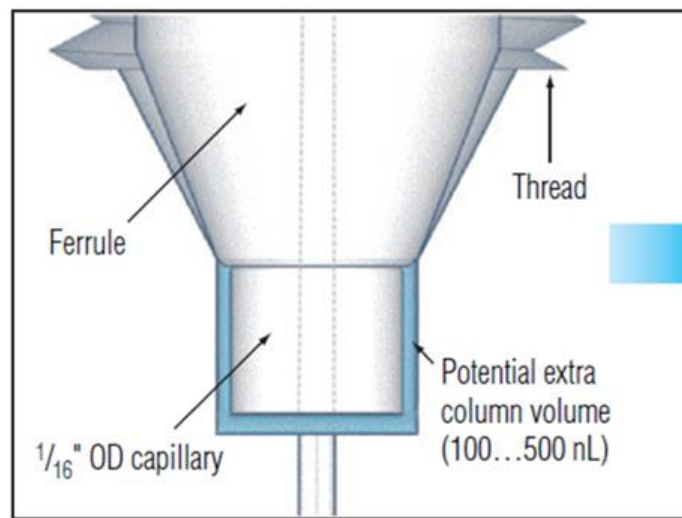
VIRTUALLY ZERO DEAD VOLUME

Seals at the tip of the tubing to minimize extra column volumes for higher efficiency and better chromatography—no more tailing peaks

UNIVERSAL CONNECTIONS

Compatible with most columns and valves and small enough to fit in the most confined spaces, such as on 10-port valves

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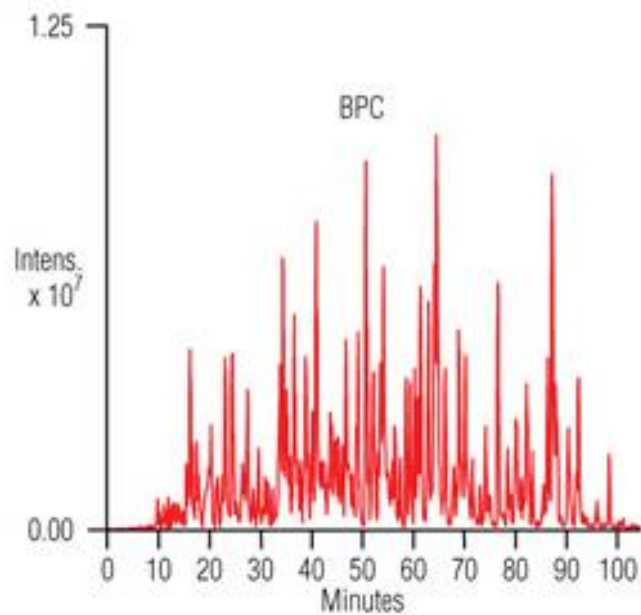
Tryptic Digest of a Protein Mixture on a Pepmap™ C18 Nanocolumn

Nanocolumn: 75 μ m i.d. x 15 cm,
PepMap100 C18, 3 μ m, 100 Å

Temperature: Ambient

Flow Rate: 300 nL/min

Solvents: (A) H₂O/acetonitrile (98:2 v/v),
0.1 % formic acid
(B) H₂O/acetonitrile (20:80 v/v),
0.08 formic acid

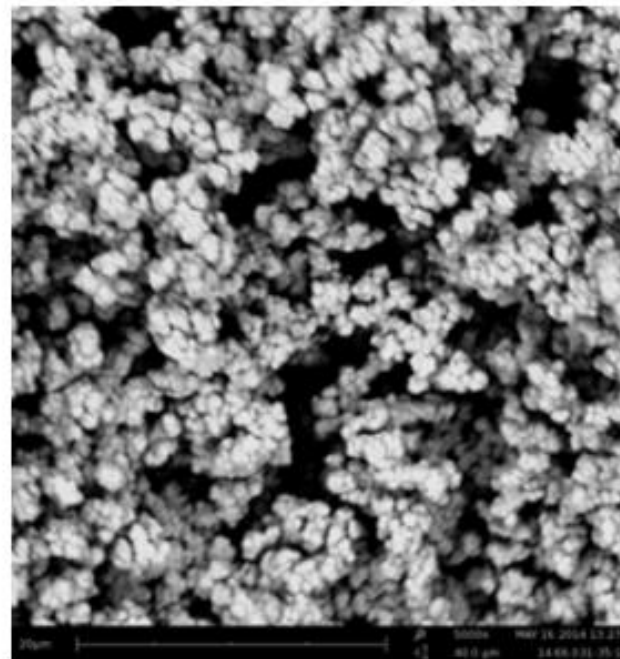
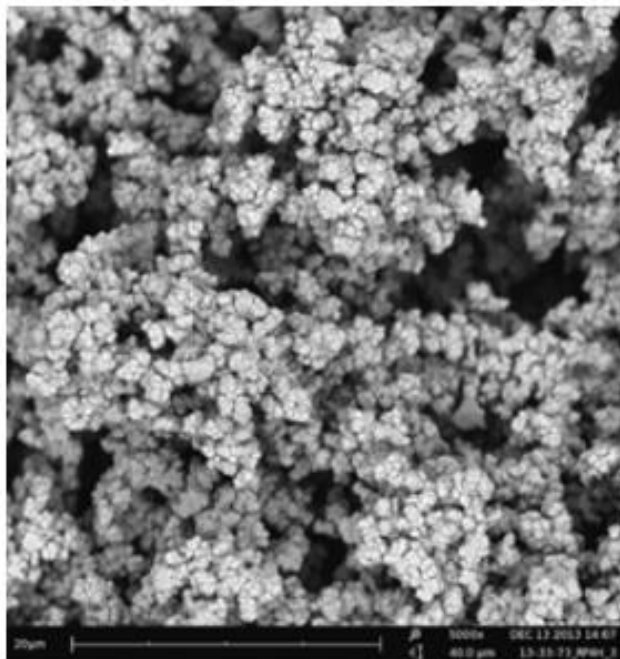
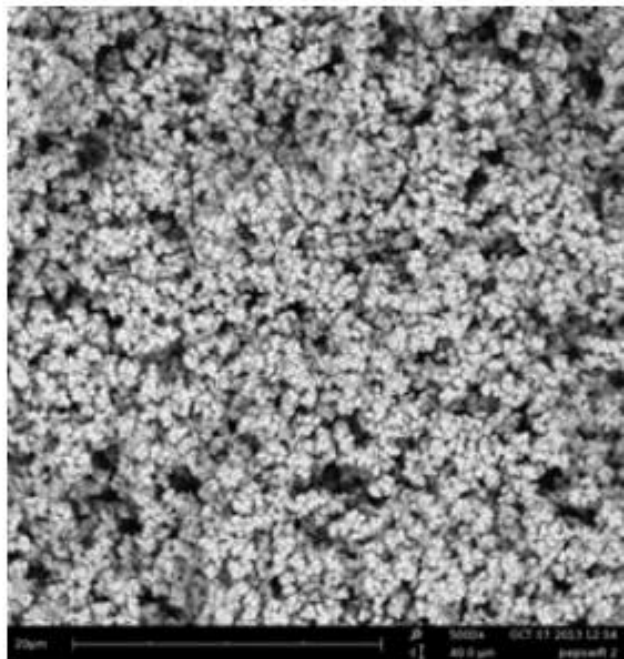
Gradient: 0–40% B in 120 min

MS: esquire3000 (Bruker)

Sample: Tryptic digest of a protein mixture:
cytochrome c, lysozyme,
alcohol dehydrogenase,
bovine serum albumin,
apotransferrin, β -galactosidase,
500 fmol each


PepSwift, ProSwift

[Více o PepSwift >>>](#)

[Více o ProSwift >>>](#)

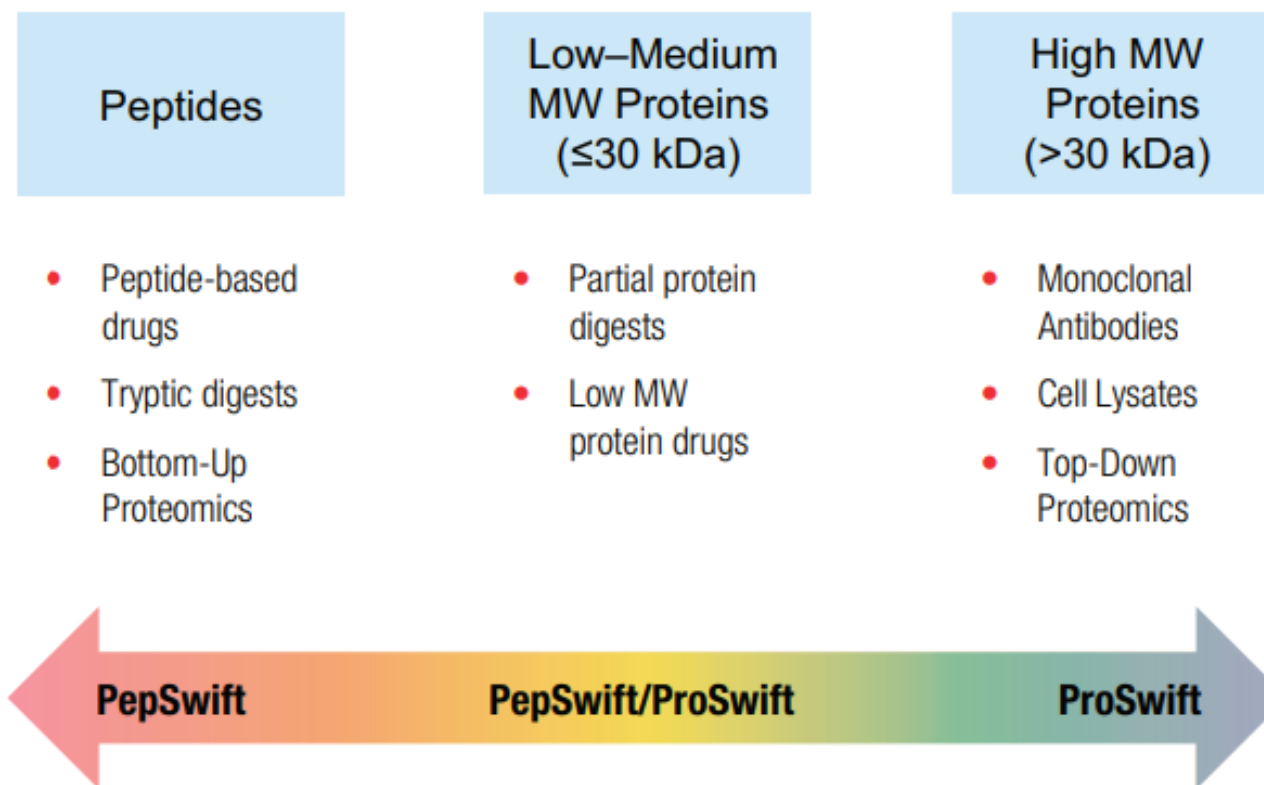
PepSwift, ProSwift

Sample: 25 µg/mL pNP-labeled sugars

1. α-glucose
2. β-glucose
3. Unknown impurity
4. α-mannose

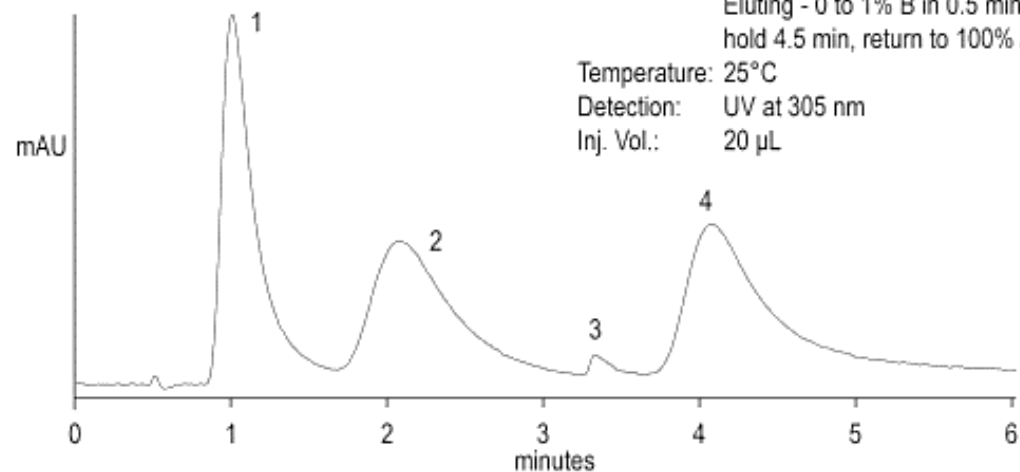Column: ProSwift ConA-1S (5 x 50 mm)
Flow Rate: 1 mL/min
Eluents: A. 50mM NaOAc, 0.2M NaCl, pH 5.3, 1mM CaCl₂
B. 100mM α-methyl mannopyranoside in eluent A
Gradient: Binding - 100% A for 1 min
Eluting - 0 to 1% B in 0.5 min, hold 4.5 min, return to 100% A
Temperature: 25°C
Detection: UV at 305 nm
Inj. Vol.: 20 µL

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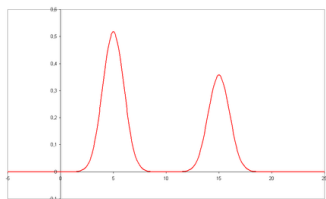
Vysoká
citlivost

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
ThermoFisher
SCIENTIFIC

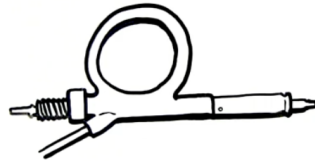thermofisher.com/lowflowHPLCcolumns

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MabPac – ProPac – PepMap – EasySpray – Acclaim

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